Operating Instructions | Betriebsanleitung | Mode d'emploi |
Instrucciones de manejo | Manuale d'uso | Handleiding

Original Operating Instructions | Original-Betriebsanleitung | Mode d'emploi original |
Instrucciones de manejo originales | Manuale d'uso originale | Originele handleiding

Flexsafe® Pro Mixer Drive Unit

Drive Unit for Use with Flexsafe® Mixing Bags and Pallettanks for Mixing

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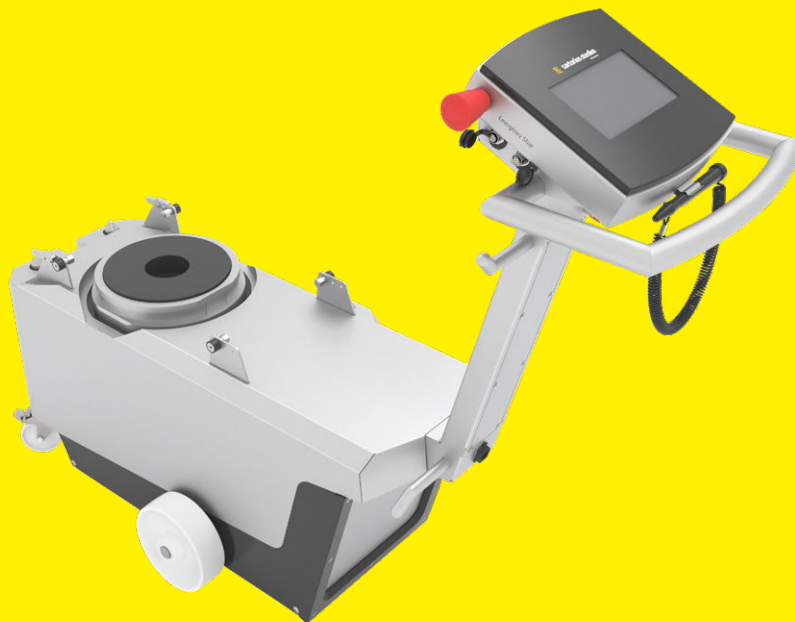
Contents

1	About these Instructions	7
1.1	Scope	7
1.2	Other Applicable Documents	7
1.3	Target Groups	7
1.4	Symbols Used	8
1.4.1	Warnings in Operating Instructions	8
1.4.2	Other Symbols	8
2	Safety Precautions	8
2.1	Intended Use	8
2.1.1	Modifications to the Device	9
2.1.2	Repairs and Maintenance Work on the Device	9
2.1.3	Foreseeable Misuse	9
2.2	Personnel Qualification	10
2.3	Significance of these Instructions	10
2.4	Functionality of the Device	10
2.5	Safety Equipment	10
2.6	Safety Information on the Device	11
2.7	Electrical Equipment	11
2.7.1	Damage to the Device's Electrical Equipment	11
2.7.2	Power Supply Cable	11
2.7.3	Potential Equalization Cable	11
2.8	Conduct in an Emergency	12
2.9	Accessories, Consumables, and Spare Parts	12
2.10	Personal Protective Equipment	12
2.11	Working on the Device	13
2.12	Order of the Required Steps	13
2.13	Moving the Device	13
2.14	Removing the Device from the Palletank® for Mixing	13
2.15	Moving Parts	14
2.15.1	Motor	14
2.15.2	Alignment Pins	14
2.15.3	Guide Rollers	14
2.16	Magnetic Forces	15
2.16.1	Effects on Implants From Magnetic Forces	15
2.16.2	Danger of Crushing Due to Magnetic Forces	15
2.17	Trip Hazard Due to Connection Cables	15
3	Device Description	16
3.1	Device Overview	16
3.2	Components for Coupling with the Palletank® for Mixing and Mixing Bag	17
3.3	Fan	17
3.4	Connections, Cable Reel and On-Off Switch	18
3.5	Coupling and Decoupling the Device on the Palletank® for Mixing	19
3.6	Acoustic Signals	19
3.7	Emergency Shut-off Switch	20
3.8	Symbols on the Device	20

4	Operating Concept	21
4.1	Display in the Main Menu	21
4.2	Display in the Mixing Process	22
4.3	Keypad	22
4.4	User Guide	23
4.5	Message	24
4.6	Recipes	24
4.7	User Administration	24
4.8	Buttons for Navigation and for Making Changes	24
4.9	Menu Structure	26
4.9.1	Menu Structure in the Program Menu	26
4.9.2	Menu Structure in the Administration Menu	26
4.10	Navigating the Menus	27
5	Installation	28
5.1	Scope of Delivery	28
5.2	Unpacking the Device	28
5.3	Prerequisites for Assembly at Installation Site	28
5.4	Removing the Transport Lock	29
6	Getting Started	30
6.1	Connecting the Power Supply	30
6.1.1	Connecting a Potential Equalization Cable	30
6.1.2	Connecting the Power Supply Cable	30
6.2	Acclimatizing the Device	31
6.3	Switching the Device On or Off	31
6.4	Moving the Device to the Installation Site	32
6.5	Connecting the Connection Cable for Remote Operation	32
7	System Settings	33
7.1	Logging Into the Device	33
7.2	Changing Regional Settings	33
7.3	Creating, Changing, or Deleting Users	34
7.3.1	Overview of User Groups	34
7.3.2	Calling up the List of Saved Users	34
7.3.3	Creating Users	35
7.3.4	Changing Users	35
7.3.5	Deleting Users	36
7.4	Modifying Access Restrictions for the Mixing Process	36
7.5	Changing the Recipe (Only "Mixing with recipe" Operating Mode)	36
7.5.1	Calling Up and Checking the Recipe	36
7.5.2	Modifying Recipe Entries	37
7.6	Parameter List	38
7.6.1	Parameters in the Program Menu	38
7.6.2	Parameters in the "Settings" "Date & time" and "Language" Menu	38
7.6.3	Parameters in the "Settings" "Access level" Menu	38

8	Operation	39
8.1	Setting Up Process Data Acquisition for USB Mass Storage	39
8.2	Selecting the Operating Mode	39
8.3	Selecting the Recipe (Only "Mixing with recipe" Operating Mode)	39
8.4	Installing the Mixing Bag in the Palletank® for Mixing	40
8.5	Inserting the Device under the Palletank® for Mixing	40
8.6	Coupling the Device to the Palletank® for Mixing and Impeller	42
8.7	Filling the Mixing Bag	43
8.8	Set the Mixing Speed and Duration (Only "Manual mixing" or "Remote mixing" Operating Mode)	43
8.9	Carrying Out the Mixing Process	44
8.9.1	Starting the Mixing Process	44
8.9.2	Pausing, Stopping, or Canceling the Mixing Process	44
8.10	Processing Messages	44
8.11	Viewing Recorded Messages	45
8.12	Ending the Mixing Process	45
8.12.1	Decoupling the Device from the Impeller and Palletank® for Mixing	45
8.12.2	Emptying and Removing the Mixing Bag	46
8.12.3	Removing the Device from the Palletank® for Mixing	46
8.13	Ending the Process Data Acquisition	47
9	Cleaning and Maintenance	48
9.1	Cleaning Device Surfaces	48
9.2	Maintenance Schedule	48
9.3	Restarting the System	48
10	Malfunctions	49
10.3	Faults when Operating the Device	52
10.4	Leaked Liquids	52
10.5	Checking the Position of the Mixing Bag and Impeller	52
10.6	Checking the Coupling with the Mixing Bag and Impeller	52
11	Decommissioning	53
11.1	Decommissioning the Device	53
12	Transport	54
12.1	Installing the Transport Lock	54
13	Storage and Shipping	55
13.1	Storage	55
13.2	Returning Devices and Parts	55
14	Disposal	56
14.1	Information on Decontamination	56
14.2	Disposing of the Device and Parts	56
14.2.1	Information on Disposal	56
14.2.2	Disposal	56

15	Technical Data	57
15.1	Dimensions and Weight	57
15.2	Ambient Conditions	57
15.3	Electrical Data	57
15.3.1	Power Supply	57
15.3.2	Electromagnetic Compatibility	58
15.4	Suitability of Palletank® for Mixing and Mixing Bags	58
15.5	Mixing Process	58
15.6	Remote Operation	59
15.7	Recipe Storage and Message Memory	59
15.8	Materials	59
15.9	Acoustic Emission	60
16	Accessories	60
17	Sartorius Service	60
18	Conformity Documents	60

1 About these Instructions

1.1 Scope

These instructions are part of the device. These instructions apply to the device in the following versions:

Device	Type
Flexsafe® Pro Mixer Drive Unit	FMD300001

1.2 Other Applicable Documents

- ▶ In addition to these instructions, observe the following documentation:
 - Palletank® for Mixing Operating Instructions
 - Information for handling mixing systems, e.g. „Flexsafe® Pro Mixer Global Operating Handling“
 - If the unit is operated remotely: Operating Instructions for Remote Control

1.3 Target Groups

These instructions are addressed to the following target groups. The target groups must possess the specified knowledge.

Target group	Knowledge and qualifications
User	The user is responsible for operating the device and for the associated work processes. They understand the hazards which may arise when working with the device and know how to prevent them. They have been trained in the operation of the device.
Operator	The operator of the device is responsible for compliance with safety requirements and workplace safety regulations. The operator must ensure that all persons who work with the device have access to the relevant information and are trained in working with the device.

1.4 Symbols Used

1.4.1 Warnings in Operating Instructions

WARNING

Denotes a danger with the risk that death or severe injury may result if it is **not** avoided.

CAUTION

Denotes a hazard that may result in moderate or minor injury if it is **not** avoided.

NOTICE

Denotes a danger with the risk that property damage may result if it is **not** avoided.

1.4.2 Other Symbols

- ▶ Required action: Describes actions which must be carried out.
- ▷ Result: Describes the result of the actions carried out.
- [] Text inside brackets refers to control and display items.
- [] Text inside brackets indicates status, warning, and error messages.

Figures on the Operating Display

The figures on the operating display of the device may deviate from those in these instructions.

2 Safety Precautions

2.1 Intended Use

The Flexsafe® Pro Mixer Drive Unit is used for controlling the impeller in a Flexsafe® Pro Mixer Bag (mixing bag). The mixing bag must be inserted into a Palletank® for Mixing.

The device must be coupled with the Palletank® for Mixing and the impeller in the mixing bag. During coupling, the device's motor surrounds the impeller in the mixing bag and drives it using magnetic force. This allows the liquids and | or solids in the mixing bag to be mixed, i.e. for processing materials.

The Palletank® for Mixing and the mixing bag must be approved for operation with the device (see Chapter "15.4 Suitability of Palletank® for Mixing and Mixing Bags", page 58).

The device can be connected to a remote control. This allows the device's individual functions to be controlled remotely, i.e. starting or pausing the mixing process. The remote control must be suitable for operation with the device (suitability, see Chapter "15.6 Remote Operation", page 59).

The device is intended exclusively for use in accordance with these instructions. Any further use beyond this is considered **improper**.

If the device is **not** used properly: The protective systems of the device may be impaired. This can lead to unforeseeable personal injury or property damage.

Operating Conditions for the Device

Do **not** use the device in potentially explosive environments. The device may only be used indoors.

The device may only be used with the equipment and under the operating conditions described in the Technical Data section of these instructions.

2.1.1 Modifications to the Device

If the device has been modified, e.g. by adding additional components: The safety of the device may be impaired or the device compliance may lose its validity.

If you have questions about modifications to the device, please contact Sartorius Stedim Biotech.

2.1.2 Repairs and Maintenance Work on the Device

Specialist knowledge about the device is required in order to carry out repair and maintenance work on it. If the device is **not** repaired or serviced by a specialist: The safety of the device may be impaired or the approval marks may lose their validity. **Do not** open the device yourself.

We recommend that any repair work, even that **not** covered by the warranty, is carried out by Sartorius Service or after consulting with Sartorius Service.

Only carry out the maintenance work described in these instructions. For maintenance work that has to be carried out by Sartorius Service, please contact Sartorius Service.

2.1.3 Foreseeable Misuse

It is only safe to use the device when operated in accordance with its intended use. The following applications, for example, are not permitted:

- Use of unsuitable or inadequately dimensioned spare parts, such as power supply cables.
- Switching on the device without the potential equalization cable having been grounded.

2.2 Personnel Qualification

If people who **do not** have sufficient knowledge on the safe handling of the device carry out work on the device: Those people may injure themselves or other people nearby.

- ▶ Ensure that all persons working on the device possess the necessary knowledge and qualifications (description see Chapter “1.3 Target Groups”, page 7).
- ▶ If a particular qualification is indicated for the actions described: Have these activities carried out by the required target group.
- ▶ If **no** particular qualification is required for the activities described: Have these activities carried out by the “user” target group.

2.3 Significance of these Instructions

Failure to follow the instructions in this manual can have serious consequences, e.g., exposure of individuals to electrical, mechanical, or chemical hazards.

- ▶ Before working with the device: Read the instructions carefully and completely.
- ▶ If these instructions are lost: Request a replacement or download the latest version from the Sartorius website (www.sartorius.com).
- ▶ Retain the instructions.
- ▶ The information contained in these instructions must be available to all individuals working with the device.

2.4 Functionality of the Device

A damaged device or worn parts may lead to malfunctions or cause hazards which are difficult to recognize.

- ▶ Only operate the device when it is safe and in perfect working order.
- ▶ Observe the maintenance intervals (intervals and maintenance work see Chapter “9.2 Maintenance Schedule”, page 48).
- ▶ Have any malfunctions or damage repaired immediately by Sartorius Service.
- ▶ Ensure the symbols that are stuck onto the device are legible.

2.5 Safety Equipment

The safety equipment on the device protects persons who work with the device against the hazards and harms associated with it, e.g., electrical current. If the device’s safety equipment is detached or modified: People may be seriously injured.

- ▶ Do **not** dismantle, or modify the safety equipment (safety equipment see Chapter “3.7 Emergency Shut-off Switch”, page 20).

2.6 Safety Information on the Device

Symbols, e.g. warnings, safety stickers, are safety information for the handling of the device. Missing safety information may lead to serious injuries.

- ▶ Do **not** remove or modify the symbols.
- ▶ Replace the symbols if they become illegible.

2.7 Electrical Equipment

2.7.1 Damage to the Device's Electrical Equipment

Damage to the device's electrical equipment, e.g. damaged insulation, can be life-threatening. There is a danger to life from contact with live parts.

- ▶ If the electrical equipment of the device is defective, cut off the power supply and contact Sartorius Service.
- ▶ Keep live parts away from moisture. Moisture can cause short-circuits.

2.7.2 Power Supply Cable

The use of an unsuitable or inadequately dimensioned power supply cable may have an unforeseeable impact and may cause life-threatening injuries as a result of electric shocks, for example.

- ▶ Only use the original power supply cable.
- ▶ If the power supply cable must be replaced: Contact Sartorius Service. Do **not** repair or modify the power supply cable.

2.7.3 Potential Equalization Cable

The lack of a potential equalization cable or the use of a defective potential equalization cable may result in damage and may cause life-threatening injuries, e.g. as a result of electric shocks.

- ▶ Only switch the device on when the potential equalization cable is connected.
- ▶ Use the potential equalization cable provided.
- ▶ If the potential equalization cable needs to be replaced: Please contact Sartorius Service. Do **not** repair or modify the potential equalization cable.

2.8 Conduct in an Emergency

If there is immediate danger of personal injury or equipment damage, e.g., due to malfunctions or dangerous situations, the device must be immediately taken out of operation.

- ▶ Press the emergency shut-off switch. This means that the power supply to the device is interrupted and all lifting or mixing motions will be stopped.
- ▶ Leave the emergency switch-off switch in the engaged position until the faults or malfunctions have been resolved.
- ▶ Malfunctions should be resolved by Sartorius Service.
- ▶ When the fault or malfunction is repaired: Disengage the emergency shut-off switch. To do so, pull out the emergency shut-off switch until the green ring on the emergency shut-off switch is visible.

2.9 Accessories, Consumables, and Spare Parts

The use of unsuitable accessories, consumables, and spare parts can affect the functionality and safety of the device and have the following consequences:

- Risk of injury to persons
 - Damage to the device
 - Malfunction of the device
 - Device failure
- ▶ You should only use approved accessories, consumables, and spare parts supplied by Sartorius Stedim Biotech.
 - ▶ Only use accessories, consumables, and spare parts that are in proper working order.

2.10 Personal Protective Equipment

Personal protective equipment protects against risks arising from the device and the materials being processed.

Protective Equipment Designation	Explanation Examples
Safety boots	Protect against injuries to the feet caused by mechanical effects.

- ▶ Wear appropriate personal protective equipment.
- ▶ When the workplace or the process, in which the device is used, requires additional safety precautions: Additionally wear the appropriate personal protective equipment.

2.11 Working on the Device

Not all areas of the device or Palletank® for Mixing can always be surveyed in full, the visibility of certain areas may be restricted, i.e. if the device is inserted or removed at the Palletank® for Mixing.

If the unit is operated while there are people in the device's operating area: People may find themselves in the danger zone between the device and the Palletank® for Mixing and are at risk of injury.

- ▶ When carrying out any work on the device, ensure beforehand that there are **no** people between the device and Palletank® for Mixing.
- ▶ Precisely define responsibilities and working areas on the device.
- ▶ Wear personal protective equipment.

2.12 Order of the Required Steps

Conducting work on the device in the incorrect order may result in unforeseen hazards. There is a risk of injury to personnel.

- ▶ When working on the device, e.g. to insert the device on the Palletank® for Mixing or for filling the mixing bag, follow the order of the work steps described in these instructions.

2.13 Moving the Device

When moving the device, there is a risk of injury, e.g. hitting someone with the device.

- ▶ When moving the device ensure that **no** personnel or objects are in the way.
- ▶ Wear personal protective equipment.

2.14 Removing the Device from the Palletank® for Mixing

If the device has been removed from the Palletank® for Mixing and comes into contact with people or people reach into the danger zone between the device and Palletank® for Mixing: There is a risk of people becoming injured, e.g. feet or hands being crushed.

- ▶ When removing the device from the Palletank® for Mixing, ensure that there are **no** people in the operating area of the device or Palletank® for Mixing.
- ▶ Do **not** reach into the area between the device and the Palletank® for Mixing while the device is being removed from the Palletank® for Mixing.
- ▶ Slowly remove the device from the guide rails of the Palletank® for Mixing. The device must be removed from the Palletank® for Mixing in a controlled manner.
- ▶ Wear personal protective equipment.

2.15 Moving Parts

2.15.1 Motor

If the user of the device or other people nearby come into contact with the motor during this process: There is a risk of fingers becoming trapped and crushed between the motor and the Palletank® for Mixing.

- ▶ Do **not** reach into the area between the device and the Palletank® for Mixing while the motor is raised.
- ▶ Ensure that there are **no** people in the operating area of the device who can reach between the motor and Palletank® for Mixing.

2.15.2 Alignment Pins

When coupling the device to the Palletank® for Mixing, the alignment pins are raised and connect to the alignment plate of the Palletank® for Mixing. If the user or other people nearby come into contact with the alignment pins or the alignment plate during this process: There is a risk of fingers becoming trapped and crushed between the alignment pins and the alignment plate.

- ▶ Do **not** reach into the area between the device and the Palletank® for Mixing while the alignment pins are raised.
- ▶ Ensure that there are **no** people in the operating area of the device who can reach between the alignment pins and alignment plate.

2.15.3 Guide Rollers

As the device is being inserted or removed, the device's guide rollers glide over the guide rails of the Palletank® for Mixing. If the user of the device or other people nearby come into contact with the guide rollers: There is a risk of fingers becoming trapped and crushed between the guide rollers and the guide rails.

- ▶ When inserting or removing the device on the Palletank® for Mixing, do **not** reach into the area between the guide rollers and guide rails of the Palletank® for Mixing.
- ▶ Ensure that there are **no** people in the operating area of the device who can reach between the guide rollers and guide rails.

2.16 Magnetic Forces

2.16.1 Effects on Implants From Magnetic Forces

The device's motor emits strong magnetic forces during operation. Magnetic forces can affect implants and other magnetic devices. People with implants can be lethally or seriously injured by these forces; magnetic or magnetized devices can also be damaged by these forces.

- ▶ Ensure that active implant carriers e.g. those with pacemakers, cochlear implants, insulin pumps, maintain a safe distance of 0.5 m from the device.
- ▶ Ensure that passive implant carriers maintain a safe distance of 0.1 m from the device.
- ▶ Magnetic or magnetized objects should **not** be stored or used near the device.

2.16.2 Danger of Crushing Due to Magnetic Forces

When the mixing bag is installed in the Palletank[®] for Mixing: High magnetic forces exist between the motor of the device and the impeller in the mixing bag. If the user of the device or other people nearby come into contact with the motor or the impeller in the mixing bag during this process: There is a risk of fingers becoming trapped and crushed.

- ▶ When working to install the mixing bag into the Palletank[®] for Mixing, observe the information in the device's operating display. The device's user guide displays the required work steps for installing the mixing bag in sequential order.
- ▶ The mixing bag must only be installed in the Palletank[®] for Mixing when the device motor is decoupled from the Palletank[®] for Mixing (the motor is fully inserted).

2.17 Trip Hazard Due to Connection Cables

If the device's connection cable, i.e. the power supply cable, is accidentally misplaced or **not** considered when moving the device: Users of the device or other people nearby might trip over the connection cable and injure themselves.

- ▶ While moving the device pay attention to where the connection cable is.

3 Device Description

3.1 Device Overview

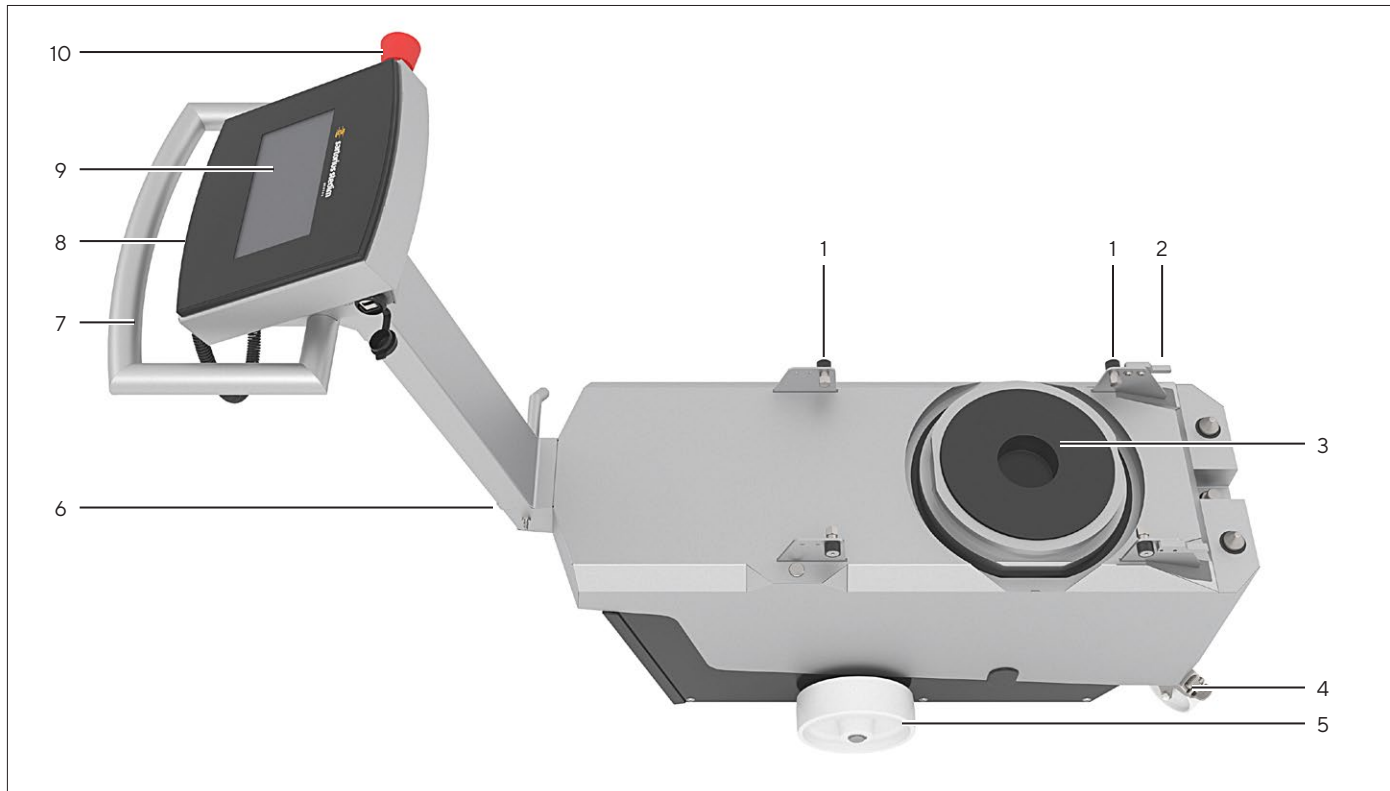
Fig.1: Flexsafe® Pro Mixer Drive unit (viewed from above)

Pos.	Designation	Description
1	Guide roller	Used to insert and remove the device underneath the Palletank® for Mixing. Glides across the Palletank® for mixing guide rails.
2	End piece	Used to ensure that the device is correctly positioned on the guide rails of the Palletank® for Mixing.
3	Motor	
4	Castors with locking brakes	Used to steer the device. Can pivot.
5	Castors without locking brakes	
6	Manufacturer's ID label	
7	Handle	Used to move and lift the device.
8	Stylus	Used to operate the operating display, e.g. when wearing gloves.
9	Operating display	Used to operate the device and control the mixing process.
10	Emergency shut-off switch	

3.2 Components for Coupling with the Palletank® for Mixing and Mixing Bag

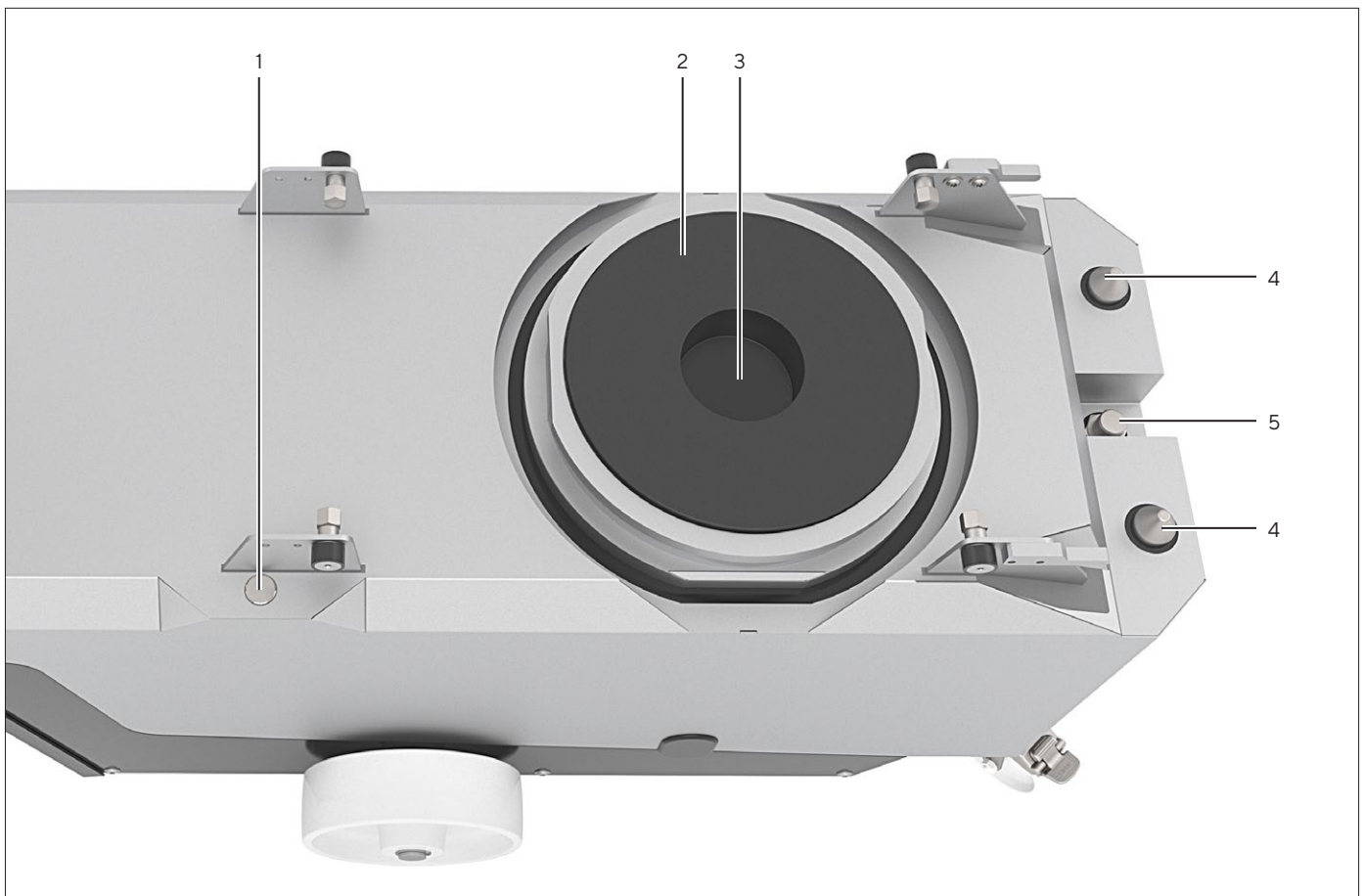
Fig.2: Components for Coupling with the Palletank® for Mixing and mixing bag (moving parts lowered)

Pos.	Designation	Description
1	Palletank® recognition sensor	Determines if the device is positioned under a Palletank® for Mixing.
2	Motor	Powers the impeller in the mixing bag. Is movable.
3	Impeller holder	Holds the impeller of the mixing bag.
4	Alignment pins	Aligns the device when rising up to the Palletank® for Mixing. Is movable.
5	Distance sensor	Determines the distance from the Palletank® for Mixing. Is movable.

3.3 Fan

The fan is used to cool the device's motor. The fan can be found on the bottom of the device.

3.4 Connections, Cable Reel and On-Off Switch

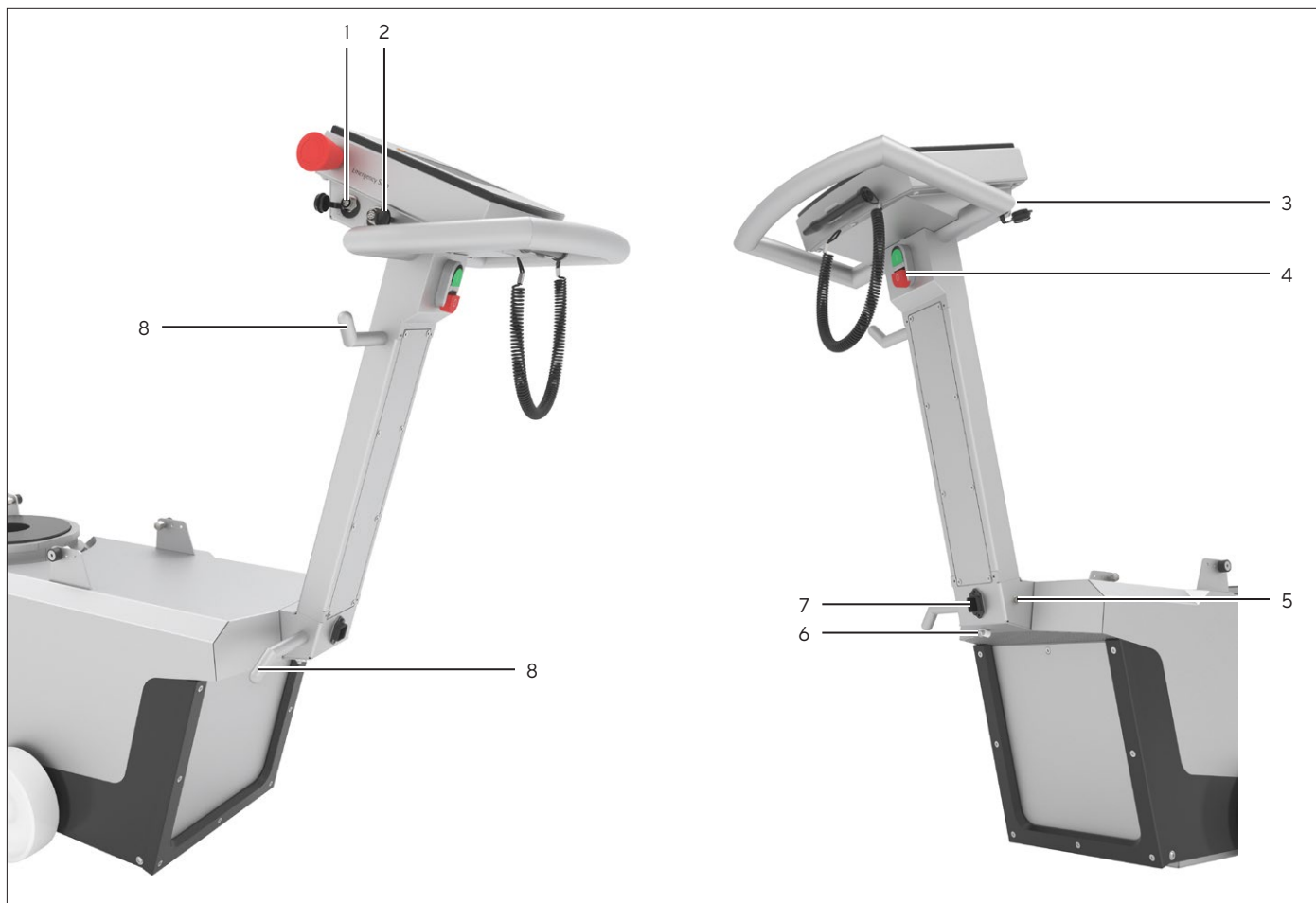
Fig. 3: Connections, cable reel and on-off switch

Pos.	Designation	Description
1	ETHERNET	Connection for Sartorius Service personnel only. Is sealed with a cover cap.
2	REMOTE	Connection for remote control. Is sealed with a cover cap.
3	USB	Connection for USB mass storage. Is sealed with a cover cap.
4	On-Off switch	Switches the device on or off.
5	Potential equalization	Connection for potential equalization cable
6	Strain relief for the power supply cable	
7	Power supply	Connection for power supply cable
8	Cable reel	Used for securing longer connection cables.

3.5 Coupling and Decoupling the Device on the Palletank® for Mixing

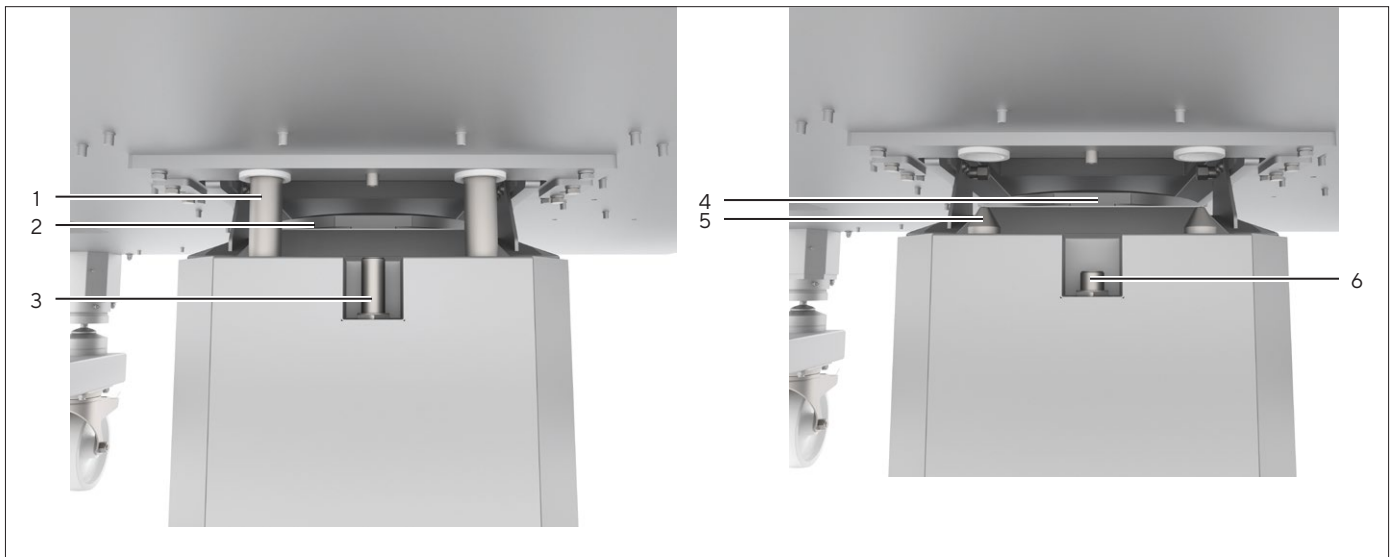
Fig. 4: Device coupled with Palletank® for Mixing and decoupled from Palletank® for Mixing (example)

Pos.	Designation	Description
1	Alignment pins coupled	The alignment pins have risen and are connected to the alignment plate of the Palletank® for Mixing.
2	Motor coupled	The motor has risen and the impeller is seated in the impeller holder of the motor.
3	Distance sensor coupled	The distance sensor has risen. The required distance from the Palletank® for Mixing has been achieved.
4	Motor decoupled	The motor has been lowered and separated from the impeller in the mixing bag.
5	Alignment pins decoupled	The alignment pins have been lowered and separated from the Palletank® for Mixing.
6	Distance sensor decoupled	The distance sensor has been lowered into the initial position.

3.6 Acoustic Signals

The device gives off acoustic signals while the motor is being raised and lowered.

3.7 Emergency Shut-off Switch

The emergency shut-off switch “EMERGENCY OFF” interrupts the device’s power supply. When the emergency shut-off switch is locked: The device **cannot** be switched on.

When the emergency shut-off switch is unlocked: The green ring on the emergency shut-off switch is visible and the device can be turned on.

3.8 Symbols on the Device

Symbol	Meaning
	During operation, parts in the device may be live. The device may only be opened by Sartorius Service. In case the electrical equipment malfunctions, the device must be immediately disconnected from the power supply.
	There is a risk of crushing when coupling the device to the Palletank® for Mixing.
	The device’s motor emits strong magnetic forces during operation.
	During operation, surfaces in the area of the motor may heat up.
	Feet may be crushed when moving the device.

4 Operating Concept

4.1 Display in the Main Menu

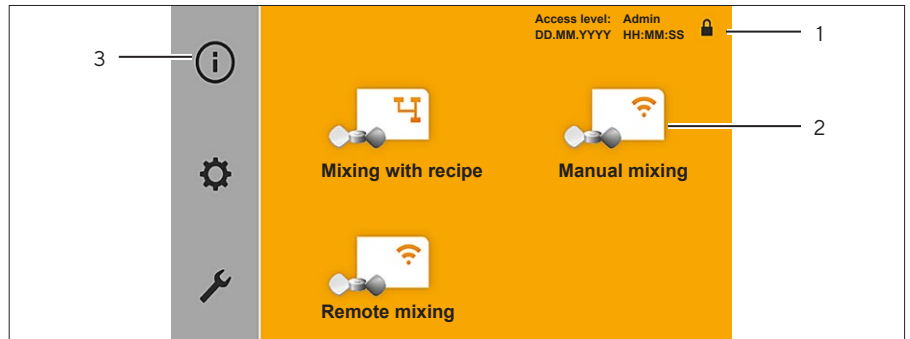
Fig. 5: "Main menu" display (example)

Pos.	Designation	Description
1	Status bar	Shows the logged-in user as well as the date and time.
2	Program menu	Used for selecting the mixing process, e.g. "Manual mixing" operating mode. The program menu display depends on the logged-in user's authorizations.
3	Administration menu	Used for performing general system settings, e.g. setting the language. The administration menu display depends on the logged-in user's authorizations.

4.2 Display in the Mixing Process

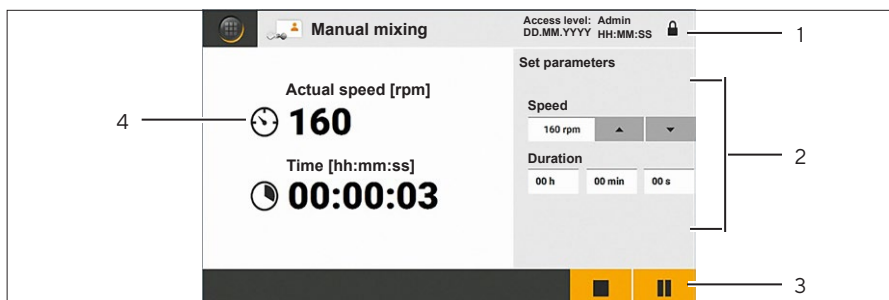
Fig. 6: Mixing process display in the “Manual mixing” operating mode (example)

Pos.	Designation	Description
1	Status bar	Displays the logged-in user, date and time, and the selected operating mode.
2	Parameter settings	Used to change the parameters for the mixing process.
3	Available functions	Shows the available functions.
4	Process data	Displays the process data for the currently running mixing process: Mixing speed and remaining or expired time.

4.3 Keypad

The keypad is used for entering values in entry fields.

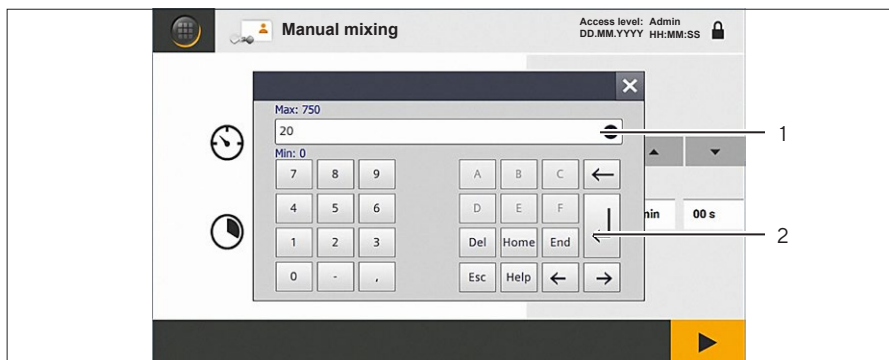
Fig. 7: Display for entering values (example)

Pos.	Designation
1	Entry field
2	Enter button

4.4 User Guide

The user guide guides the user through the process step routine and provides support with entering values in entry fields.

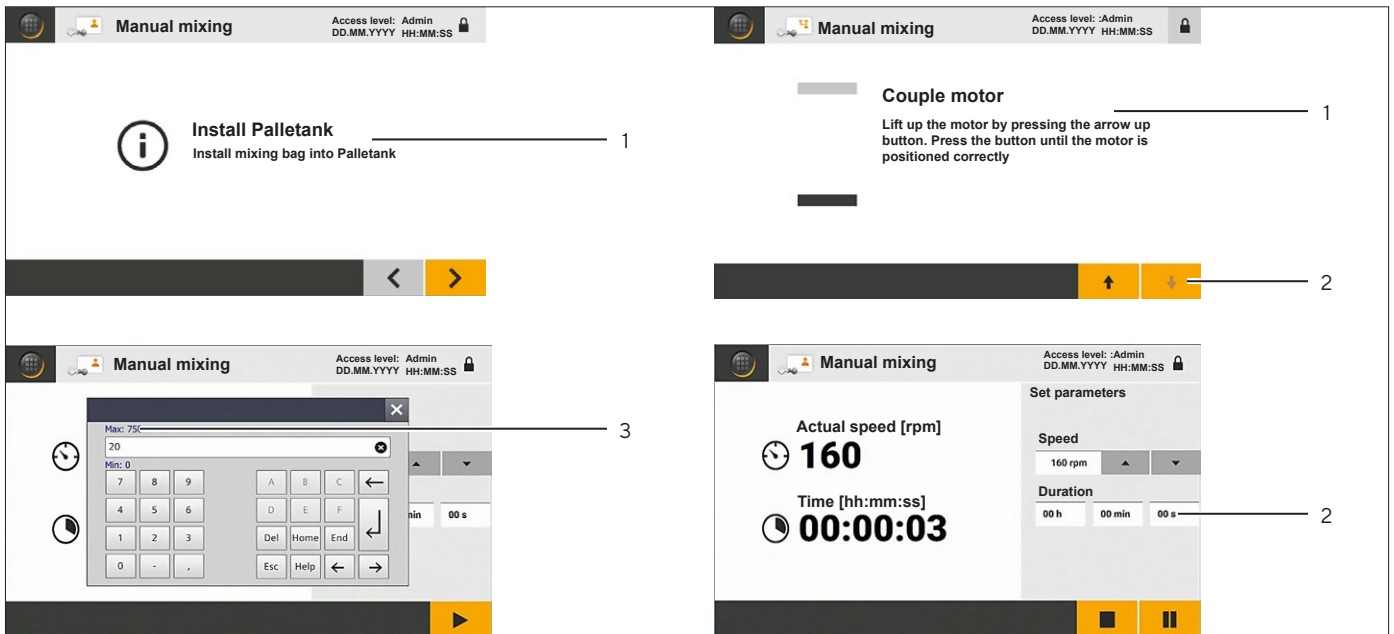
Fig. 8: User guide (example)

Pos.	Designation	Description
1	Work step	Displays the current work step in a process step routine, e.g. for coupling the motor.
2	Input assistance	Functions that are not available are grayed out, e.g. buttons or entry fields.
3	Parameter limit value	

4.5 Message

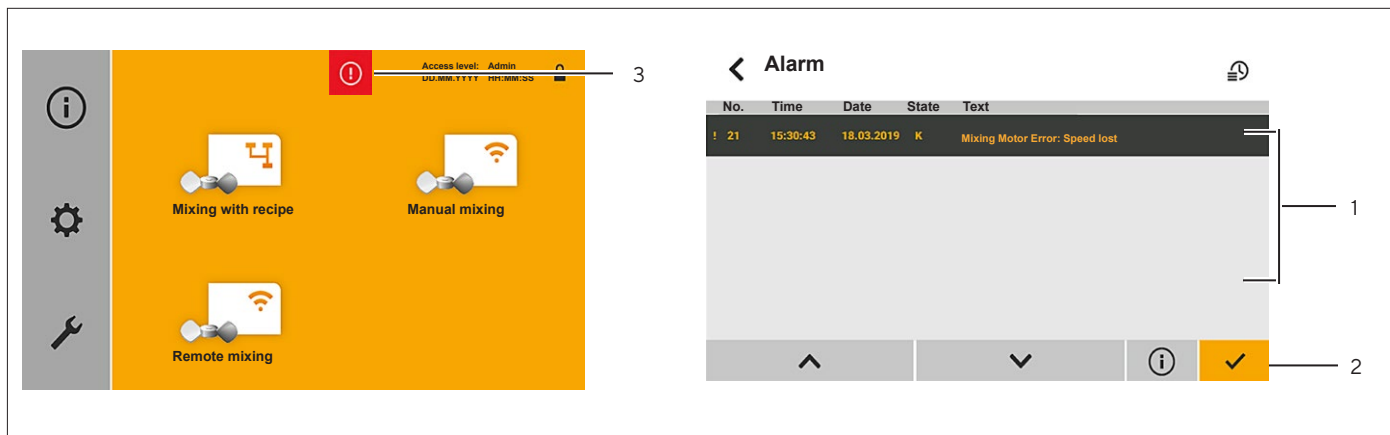
Fig. 9: Message (error message, example)

Pos.	Designation	Description
1	Error list	Displays the saved error messages and warning messages.
2	Available functions	Shows the available functions.
3	Message type	<ul style="list-style-type: none"> - Indicates that a message is present. - Defines the type of message, e.g. an error message.

4.6 Recipes

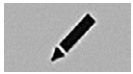
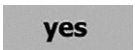
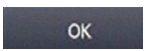
Recipes are process steps for a mixing process, which are automatically carried out. Recipes contain built-in phases for which different mixing speeds and durations can be set.

4.7 User Administration

Some functions of the device are access restricted. It is possible to create users with different authorizations.

4.8 Buttons for Navigation and for Making Changes

Pos.	Symbol	Designation	Description
1		[Home] button	Opens the main menu.
2		[Logout] button	<ul style="list-style-type: none"> - Indicates that a user is logged in on the device. - Logs the user out.
3		[Login] button	<ul style="list-style-type: none"> - Indicates that no user is logged in to the device. - Opens a dialog for logging into the device.
4		[Raise] button [Lower] button	Allows the motor and the associated moving components of the device to be raised (coupling) or lowered (decoupling).

Pos.	Symbol	Designation	Description
5		[Start] button	Starts a mixing process.
6		[Pause] button	Interrupts a mixing process.
7		[Stop] button	Stops a mixing process or the remote operation of the device.
8		[Previous display] button [Next display] button	Switches to the previous or next display in a process step routine or in a menu.
9		[Previous recipe] button [Next recipe] button	Switches to the previous or next recipe in a list of saved recipes.
10		[Value higher] button [Value lower] button	Increases or decreases the mixing speed in increments of 5.
11		[List up] button [List down] button	Scrolls up or down in a list.
12		[Details] button	Opens the detailed display of an error message or warning message.
13		[Info] button	Opens the information for an error message or warning message.
14		[Confirm] button	Confirms the settings and information in a display.
15		[Edit] button	Opens a display for editing a recipe.
16		[Error] button	<ul style="list-style-type: none"> - Indicates that an error is present. - Opens the error list.
17		[Warning] button	<ul style="list-style-type: none"> - Indicates that a warning is present. - Opens the error list.
18		[YES] button	Confirms the execution of an action in a dialog.
19		[NO] button	Cancels the execution of an action in a dialog.
20		[Close] button	Closes a display in a process step routine, a dialog, or the keypad.
21		[Cancel] button	Interrupts the entry in a dialog.
22		[OK] button	Confirms the entry in a dialog.

4.9 Menu Structure

4.9.1 Menu Structure in the Program Menu

► Navigating in menus (see Chapter “4.10 Navigating the Menus”, page 27).


Level 1	Description reference
Mixing with recipe 	The parameters for the mixing process are saved in recipes. Recipes must be called up on the device and started.
Manual mixing 	The parameters for the mixing process must be entered manually on the device.
Remote mixing 	The mixing process must not be set up and started on the device, but rather on an external control unit.

4.9.2 Menu Structure in the Administration Menu


► Navigating in menus (see Chapter “4.10 Navigating the Menus”, page 27).

Level 1	Level 2	Level 3	Description reference
System information 	Operating data		Displays recorded operating times since the device was started-up: <ul style="list-style-type: none"> - Operating time for lifting processes in minutes - Operating time for mixing processes in hours - Total operating time in hours
	Error history		Displays the error list.
	Software		Displays information on the software version of the device.
Settings 	Date & time		Set the date and time.
	Language		Set the language.
	User administration	User login <hr/> User administration	Define an access restriction for the program menu. <hr/> View and manage saved users.
Maintenance			Access only for Sartorius Service 

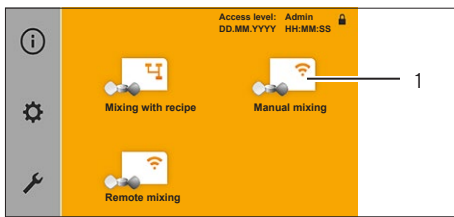
4.10 Navigating the Menus

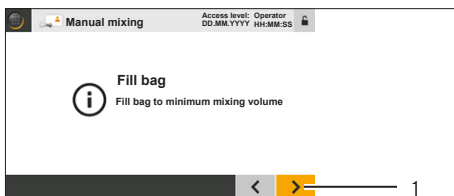
- ▶ To access a menu: Press the button for the menu, e.g. the [Manual mixing] button (1).
- ▶ The first display in the process step routine or the submenu is shown.

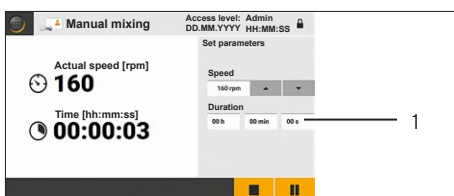- ▶ To go to the next display in a process step routine: Press the [Next display] (1) button.
- ▶ The next display for the process step routine appears.

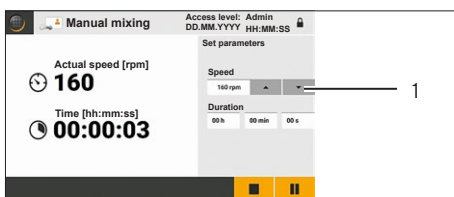- ▶ If a value needs to be entered in an entry field or changed:
 - ▶ Press the entry field (1).
 - ▶ The keypad appears.
 - ▶ Enter the desired value using the keypad.
 - ▶ Press the [Enter] button.
 - ▶ The keypad closes.
 - ▶ The entered value is displayed in the entry field.

- ▶ If a value needs to be entered in an entry field via the interface elements, e.g. the mixing speed: Hold down the button until the desired value is shown, e.g. hold down the [Value higher] or [Value lower] button until the desired value is displayed.

- ▶ To cancel a process and return to the main menu: Press the [Home] or [Close] button.

- ▶ To log out of the device: Press the [Logout] button in the status bar.
- ▶ The [Logout] button is hidden.

5 Installation

5.1 Scope of Delivery

Item	Quantity
Flexsafe® Pro Mixer Drive Unit	1
Screw to replace the transport lock screw	2
Country-specific power supply cable (EU)	1
Country-specific power supply cable (USA)	1
Potential equalization cable, combined with power supply cable and additional power supply plug	1
Operating instructions	1
Technical documentation	1

5.2 Unpacking the Device

Procedure

- ▶ Unpack the device.
- ▶ If the device is stored temporarily: Observe the storage information (see Chapter "15.2 Ambient Conditions", page 57).

5.3 Prerequisites for Assembly at Installation Site

Procedure

- ▶ Make sure that the following conditions are met at the installation site:


Condition	Features
Ambient conditions	Suitability tested (see Chapter "15.2 Ambient Conditions", page 57)
Setup surface	<ul style="list-style-type: none"> - Stable and level surface - Sufficient space for the device and for moving the device (see Chapter "15.1 Dimensions and Weight", page 57 for required dimensions) - Sufficient load-bearing capacity for the device and the peripheral devices even when full (device weight see Chapter "15.1 Dimensions and Weight", page 57; weight of the peripheral devices see instructions for the peripheral devices, e.g. Palletank® for Mixing)
Suitability of Palletank® for Mixing and mixing bag	The suitability of the Palletank® for Mixing and mixing bags has been checked (suitability, see Chapter "6.1.2 Connecting the Power Supply Cable", page 30).

5.4 Removing the Transport Lock


The device is protected with a transport lock on delivery. The transport lock consists of 2 transport lock screws. The transport lock screws must be removed in order to operate the device.

Personnel: 2 persons, user

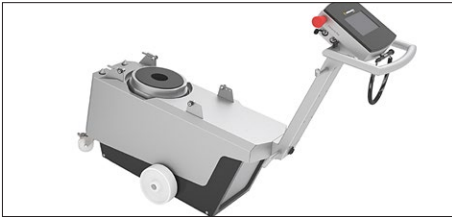
Procedure

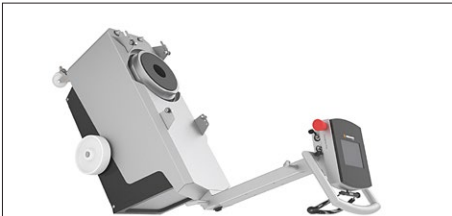- ▶ Hold the device firmly by the handle and push it down in a controlled manner.

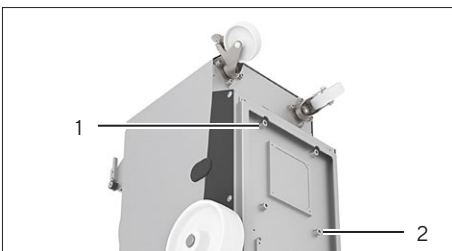- ▶ Push the device down until the handle is resting on the floor.
- ▶ Have a second person hold the device firmly by the handle, so the device remains in this position during installation

- ▶ Unscrew the transport lock screws (1) and (2). The transport lock screws are marked in red.
- ▶ Put the transport lock screws aside should the device ever need to be transported again later.
- ▶ Screw the 2 supplied screws into the device as a replacement for the removed transport lock screws.
- ▶ While holding the device by the handle, tip it downwards once again in a controlled manner so that it stands on the castors.
- ▶ **⚠ CAUTION** Risk of injury due to uncontrolled rolling away of the device! Place it on a level surface and secure the brakes. Wear personal protective equipment.

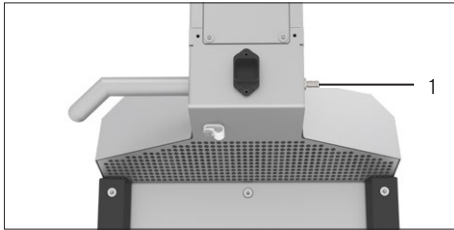
6 Getting Started

6.1 Connecting the Power Supply

6.1.1 Connecting a Potential Equalization Cable

Procedure


- ▶ **⚠ WARNING** Risk of injury due to electric shock when using an unsuitable potential equalization cable! Use the potential equalization cable provided. The potential equalization cable is combined with the power supply cable.
- ▶ Check whether the mains plug of the potential equalization cable matches the power connections at the installation site.
 - ▶ If necessary: Contact Sartorius Service.
- ▶ Connect the potential equalization cable to the “Potential equalization” connection (1).
- ▶ Connect the mains plug of the potential equalization cable to the power supply provided at the installation site.

6.1.2 Connecting the Power Supply Cable

Procedure

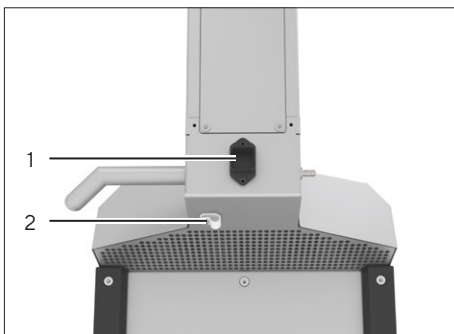- ▶ **⚠ WARNING** Risk of injury due to electric shock when using an unsuitable power supply cable! Use the power supply cable provided.
- ▶ Check whether the country-specific power plug matches the power connections at the installation site.
 - ▶ If required: Contact Sartorius Service (for available power supply cables, see Chapter “16 Accessories”, page 60).
- ▶ Unscrew the nut for the strain relief (2) and remove.
- ▶ Insert the power supply cable into the cable holder.
- ▶ Reinsert the nut for the strain relief and tighten by hand.
- ▶ Press the grooved surface on the plug of the power supply cable and connect the plug to the “power supply” connection (1).
- ▶ Connect the mains plug of the power supply cable to the power supply provided at the installation site.

6.2 Acclimatizing the Device

If the device is moved from one location to another location and the temperature at the current location is lower than the temperature at the next location: The temperature difference can lead to condensation from humidity in the device (moisture formation). Moisture in the device can lead to malfunctions. To avoid malfunctions the device must be acclimatized.

Vorgehen

- ▶ If the device was stored or used at a lower temperature than the temperature at the installation site where the device is intended to be used next: Acclimatize the device. To do so, carry out the following acclimatization steps:

Acclimatization step	Chapter, page
1. Switch on the device at the current location.	6.3, 31
2. Leave the device switched on for at least one hour.	
3. Switch the device off.	
4. Disconnect the power supply cable and potential equalization cable from the power supply at the current location and secure to the device's cable reel.	
5. Immediately move the device to the next location where the device is intended to be used.	6.4, 32
6. Immediately connect the mains plug of the power supply cable and potential equalization cable to the power supply at the installation site.	
7. Switch on the device on.	6.3, 31
8. Leave the device switched on for at least one hour. Once this waiting time has elapsed, processes can be carried out on the device.	

6.3 Switching the Device On or Off

Procedure

- ▶ **⚠ WARNING** Risk of injury due to electric shock in the event of potential equalization failure! The potential equalization cable must be connected (see Chapter "6.1.1", Page30).
- ▶ To switch the device on:
 - ▶ Press the on-off switch so that it is in the "1" position.
 - ▶ Wait until the main menu is shown.
- ▶ To switch the device off: Press the on-off switch so that it is in the "0" position.

6.4 Moving the Device to the Installation Site

Procedure

- ▶ Release the 2 locking brakes on the castors.
- ▶ **⚠ WARNING** Risk of injury while moving the device! When moving the device ensure that **no** personnel or objects are in the way. Wear personal protective equipment.
- ▶ **⚠ CAUTION** The connection cables pose a trip hazard! While moving the device pay attention to where the connecting cable is.
- ▶ Roll the device to the intended installation site.
- ▶ **⚠ CAUTION** Risk of injury due to uncontrolled rolling away of the device! Place it on a level surface and secure the brakes. Wear personal protective equipment.

6.5 Connecting the Connection Cable for Remote Operation

If the device is to be used in remote operation: A connection cable for remote operation must be connected.

Procedure

- ▶ Check whether the remote control and the connection cable for the remote control are suitable for operation of the device (see Chapter "15.6 Remote Operation", page 59).
- ▶ Connect the connection cable for remote operation to the "REMOTE" connection.
- ▶ Connect the other end of the remote operation connection cable to the external control (connection, see guide for remote control).

7 System Settings

7.1 Logging Into the Device

The device comes with preinstalled users. The user classed as “supervisor” must initially log into the device. The user classed as “supervisor” is automatically selected upon delivery. The standard password for this user is “1111”.

Procedure

- ▶ Press the [Login] button in the status bar.
- ▷ The login dialog appears.
- ▶ Enter the personal password or the standard password in the entry field.
- ▶ To confirm the entry: Press the [OK] button.
- ▷ The main menu appears.

7.2 Changing Regional Settings

Requirements

The personal user profile has the user group “group” and the password is recognized.

Procedure

- ▶ Call up the “Settings” menu in the administration menu.
- ▶ If the time setting or the date for the device needs to be changed:
 - ▶ Press the [Date & time] button.
 - ▷ The date and the time are displayed.
 - ▶ If necessary: Change the value in the entry field.
- ▶ If the language needs to be changed:
 - ▶ Press the [Language] button.
 - ▷ The currently selected language is marked with a check symbol.
 - ▶ Press the button for the desired language.

7.3 Creating, Changing, or Deleting Users

7.3.1 Overview of User Groups

Users can be created on the device. The users must be assigned to a user group. Within the user group, permissions to carry out certain functions have been defined.

To access the access restricted functions on the device: The users must log-in.

User group	Authorizations			
	View system information in the "System information" menu	Carry out mixing process (Mixing)	Change system settings (Settings)	Carry out maintenance work (Maintenance)
Users	0 or 1	0 or 1	-	-
Group	0 or 1	0 or 1	2	-
Administrator group	0 or 1	0 or 1	2	Only Sartorius Service

0 = no access restriction

1 = can be access restricted; if access restricted: Only accessible by user group "Users", "Group" or "Administrator group"

2 = is access restricted, accessible only by user group "Group" or "Administrator group"

7.3.2 Calling up the List of Saved Users

Requirements

The personal user profile has the user group "group" and the password is recognized.

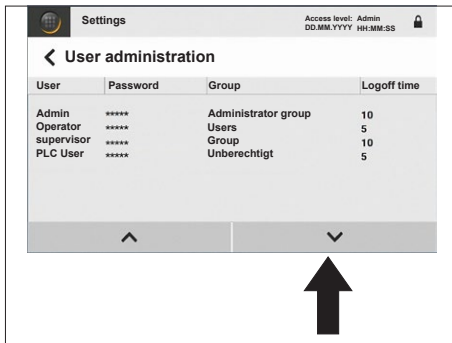
Procedure

- ▶ Call up the "Settings" menu in the administration menu.
- ▶ Press the [User administration] button.
- ▶ Press the [User administration] button.
- ▷ A list of the saved users is displayed.

7.3.3 Creating Users

Procedure

- ▶ In the list of saved users, hold down the [List down] button until the end of the list is reached.
- ▶ Double click in the empty row.
- ▷ A new entry is created.



7.3.4 Changing Users

The users classed as “supervisor” and “Operator” come pre-installed with standard passwords. We recommend that you change these standard passwords:

- Standard supervisor password: 1111
- Standard operator password: 0000

Procedure

- ▶ If a saved user needs to be changed, e.g. the password for the user:
 - ▶ Click on the corresponding user entry, e.g. on the [Password] table cell.
 - ▷ The table cell is selected.
 - ▶ Press the entry again.
 - ▷ The keypad appears.
 - ▶ Change the value in the entry field (for setting options, see Chapter “7.6.3 Parameters in the “Settings” | “Access level” Menu”, page 38).

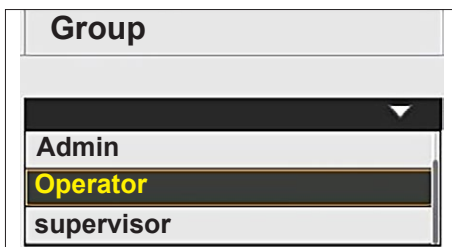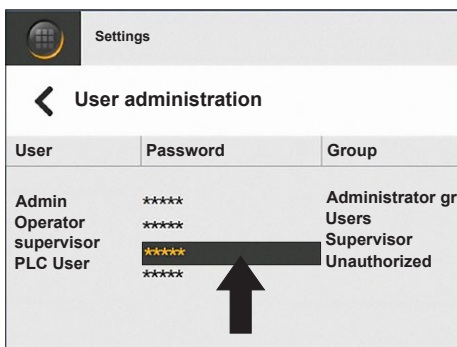- ▶ If the user group for an entry needs to be changed:
 - ▶ Press the [Group] table cell for the user.
 - ▶ The [Group] table cell is marked and an arrow symbol appears in the table cell.
 - ▶ Click on the arrow symbol.
 - ▷ A drop-down list is opened.
 - ▶ Select the desired user group. To do this, click on the user group button, e.g. the [Operator] button.
 - ▶ To close the drop-down list: Click on the arrow symbol.

7.3.5 Deleting Users

Procedure

- ▶ Go to the desired entry in list of saved users. To do so, press the [List up] or [List down] buttons.
- ▶ Double-click on the desired entry in the [User] table cell.
- ▷ The table cell is selected.
- ▶ Press the entry again.
- ▷ The keypad appears.
- ▶ Delete the entered name so that the keypad entry field is empty.
- ▶ Press the [Enter] button.
- ▷ The user is deleted from the list of saved users.

7.4 Modifying Access Restrictions for the Mixing Process

Procedure

- ▶ Call up the “Settings” menu in the administration menu.
- ▶ Press the [Access level] button.
- ▶ Press the [User login] button.
- ▷ The current access rights are marked with a check symbol.
- ▶ Change the value in the entry field (for setting options, see Chapter “7.6.3 Parameters in the “Settings” | “Access level” Menu”, page 38).

7.5 Changing the Recipe (Only “Mixing with recipe” Operating Mode)

7.5.1 Calling Up and Checking the Recipe

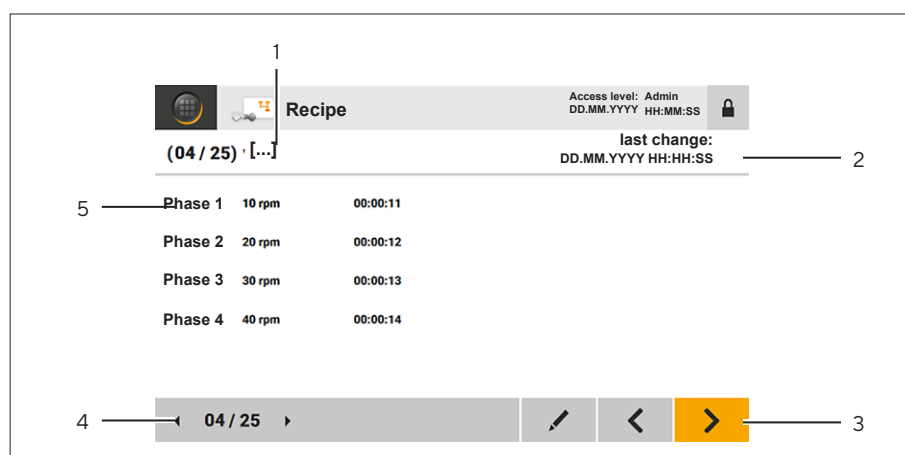
Fig.10: [Recipe] display (example)

Pos.	Designation
1	Recipe name
2	Last change (time stamp)
3	Available functions
4	List of recipes
5	Saved phase

Requirements

If the operating modes are access restricted for the mixing process: The personal user profile has the user group "Users" user group, the personal password is known.

Procedure

- ▶ Press the [Mixing with recipe] button in the program menu.
- ▷ The [Recipe] display opens.
- ▶ In the list of recipes, hold down the [Previous recipe] or [Next recipe] button until the desired recipe is displayed.

7.5.2 Modifying Recipe Entries

Procedure

- ▶ Call up the desired recipe.
- ▶ Click on the [Edit] button in the list of available functions.
- ▶ The [Edit recipe] display appears.
- ▶ Change the value in the entry fields for the recipe (for setting options, see Chapter "7.6.1 Parameters in the Program Menu", page 38).
- ▶ To save the changes: Press the [Confirm] button.
- ▶ The dialog [Are you sure you want to save?] appears.
- ▶ To save the changes to the recipe: Press the [YES] button.
- ▷ The time stamp for the recipe is updated.

7.6 Parameter List

7.6.1 Parameters in the Program Menu

Parameters	Setting values	Explanation
Speed [rpm]	20 – 750	Define the mixing speed.
Duration [00 h/00 min/00 s]	00:00:01 – 99:59:59	<ul style="list-style-type: none"> – Define the duration for a mixing process or for a phase in a recipe. – If no duration is defined: The elapsed time will be displayed in the mixing process display.
Number of phases per recipe	10	Only in the "Mixing with recipe" operating mode: <ul style="list-style-type: none"> – Define values for phases for a mixing process. – If no values are entered in a phase: The phase is not displayed in the recipe.

7.6.2 Parameters in the "Settings" | "Date & time" and "Language" Menu

Parameters	Setting values	Explanation
Date & time	DD.MM.YYYY	
	HH:MM:SS	
Language	English*	
	French	
	Spanish	
	Italian	
	German	

* Factory setting

7.6.3 Parameters in the "Settings" | "Access level" Menu

Parameters	Setting values	Explanation
User login	Yes*	Activate access restriction for the program menu.
	No	Deactivate access restriction for the program menu.
User administration		
User	User name	
Password	Password	
Group	<ul style="list-style-type: none"> – Users – Group – Administrator group 	Assign user to a user group.
Logoff time	00 – 59	Define automatic timeout value for users.

* Factory setting

8 Operation

8.1 Setting Up Process Data Acquisition for USB Mass Storage

The process data from the mixing process can be saved on a USB mass storage device (USB stick). The process data are saved in a text file. The text file can be exported and statistically evaluated, e.g. in Excel.

The following process data are included in the text file:

- Error messages or warning messages
- Speed (recorded in 1 minute intervals)
- Start and end of the mixing process
- Maximum stored values: 10.000

Procedure

- ▶ Connect a USB stick to the “USB” port.
- ▷ The process data is recorded as soon as a mixing process is started.

8.2 Selecting the Operating Mode

Requirements

If the program menu is access-restricted: The personal user profile has the user group “Users” user group, the personal password is known.

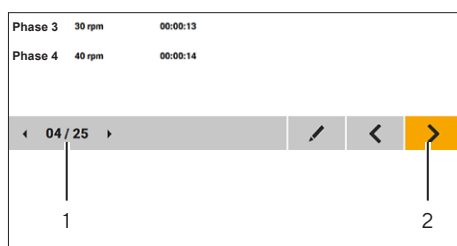
Procedure

- ▶ Switch on the device.
- ▶ Select the desired operating mode in the program menu. To do this, press the button of the operating mode, e.g. the [Mixing with recipe] button.
- ▶ If the login dialog is shown: Log-in to the device.
- ▷ The process step routine is launched. The necessary working steps are displayed one after the other.

8.3 Selecting the Recipe (Only “Mixing with recipe” Operating Mode)

Procedure

- ▶ In the list of recipes (1), hold down the [Previous recipe] or [Next recipe] button until the desired recipe is displayed.
- ▶ Check the recipe settings.
 - ▶ If necessary: Adjust the recipe, (see Chapter “8.9 Carrying Out the Mixing Process”, page 44).
- ▶ To move onwards in the process step routine: Press the [Next display] button (2).

8.4 Installing the Mixing Bag in the Palletank® for Mixing

Requirements

The [Install Palletank] display is shown in the operating display.

WARNING

Danger of crushing due to magnetic forces!

High magnetic forces exist between the motor of the device and the impeller in the mixing bag.

- ▶ Observe the information in the operating display of the device. The device's user guide displays the required work steps for installing the mixing bag and the impeller one after the other.
 - ▶ The mixing bag must only be installed in the Palletank® for Mixing when the device motor is decoupled from the Palletank® for Mixing (motor inserted).
-

Procedure

- ▶ Insert the mixing bag into the Palletank® for Mixing (see instructions for the Palletank® for Mixing).
- ▶ Confirm the process. To do so, press the [Next display] button.
- ▶ The [Magnetic shield] display appears in the user guide.
- ▶ On the underside of the Palletank® for Mixing, remove the magnetic protective cap from the impeller.
- ▶ Confirm the process. To do so, press the [Next display] button.

8.5 Inserting the Device under the Palletank® for Mixing

Requirements

- The Palletank® for Mixing is ready for operation.
- The [Dock Palletank] display is shown in the operating display.

CAUTION

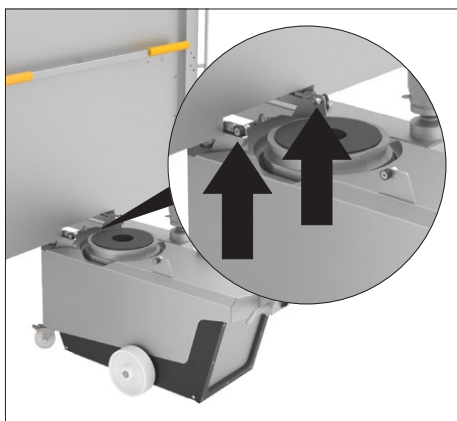
Risk of crushing due to moving parts!

When raising the device under the Palletank® for Mixing, the device's guide rollers glide over the guide rails of the Palletank® for Mixing.


- ▶ Do **not** reach into the area between the guide rollers and guide rails of the Palletank® for Mixing.
 - ▶ Ensure that there are **no** people in the operating area of the device who can reach between the guide rollers and guide rails.
-

Procedure

- ▶ Release the 2 locking brakes on the castors.
- ▶ **⚠ WARNING** Risk of injury while moving the device! When moving the device, ensure that **no** personnel or objects are in the way. Wear personal protective equipment.
- ▶ Align the device in front of the Palletank[®] for Mixing. The guide rollers must be inside of the guide rails of the Palletank[®] for Mixing.

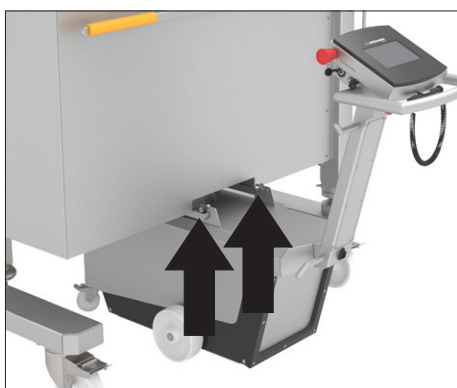- ▶ Press the handle downwards slightly so that the front guide rollers glide onto the guide rails of the Palletank[®] for Mixing.
- ▶ Push the device forwards until the rear guide rollers are in front of the guide rails of the Palletank[®] for Mixing.

- ▶ Pull the handle upwards so that the rear guide rollers glide onto the guide rails of the Palletank[®] for Mixing.

- ▶ Push the device completely under the Palletank[®] for Mixing. The device must glide over the 2 rolling elements on the guide rails of the Palletank[®] for Mixing.
- ▶ The device has been inserted under the Palletank[®] for Mixing.
- ▶ To confirm the process: Press the [Confirm] button.

8.6 Coupling the Device to the Palletank® for Mixing and Impeller

To couple the device to the Palletank® for mixing and the impeller: The motor, the alignment pins and the distance sensor must be raised.

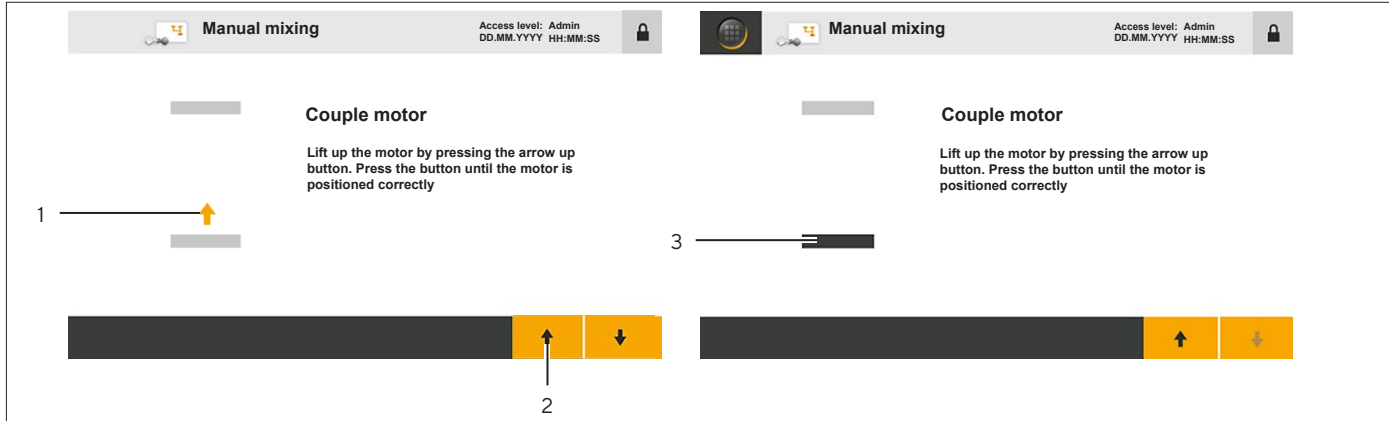
Fig. 11: Display [Couple motor] (motor in lower and upper position, example)

Pos.	Designation	Description
1	Direction of movement	Displays the direction of movement of the motor.
2	[Raise] button	If the [Raise] button is held down: The motor and the associated components are raised.
3	Positioning bar	Indicates the position of the motor: <ul style="list-style-type: none"> - Positioning bar colored black: Current position of the motor. - Positioning bar up: The motor is being raised. - Positioning bar in the middle: The motor is currently being raised or lowered. - Positioning bar down: The motor is being lowered.

Requirements

The “Couple motor” display appears in the user guide.

⚠ WARNING

Risk of crushing due to moving parts!

When coupling the device, the moving parts rise.

- ▶ Do **not** reach into the area between the device and the Palletank® for Mixing while the moving parts are rising.
- ▶ Ensure that there are **no** people in the operating area of the device who can reach between the moving parts and Palletank® for Mixing.

Procedure

- ▶ Check the position of the motor in the “Couple motor” display. The positioning bars must be in the lowest position.
- ▶ If the positioning bars are **not** in the lowest position: Click on the [Lower] button and hold it down. This causes the motor and the associated components to be lowered to the starting position.
- ▶ Click on the [Raise] button and hold it down.
- ▶ The device connects itself with the Palletank® for Mixing and the impeller in the mixing bag. During this process the distance sensor, the alignment pins, and the motor rise and the device emits acoustic signals.

8.7 Filling the Mixing Bag

Requirements

The [Fill bag] display is shown in the operating display.

Procedure

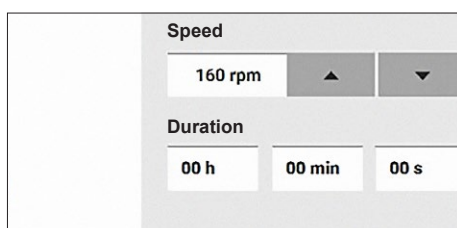
- ▶ Fill the mixing bag (see instructions for the Palletank® for Mixing).
- ▶ To confirm the process: Press the [Next display] button.

8.8 Set the Mixing Speed and Duration (Only “Manual mixing” or “Remote mixing” Operating Mode)

Requirements

The user guide shows the “mixing process” display.

Procedure

- ▶ If the device is operated in the “Manual mixing” operating mode: Specify the desired value for the mixing speed and duration (for setting options, see Chapter “7.6.1 Parameters in the Program Menu”, page 38).
- ▶ If the device is operated in the “Remote mixing” operating mode: Set the desired value for the mixing speed and duration on the remote control unit (see the remote control unit instructions).

8.9 Carrying Out the Mixing Process

8.9.1 Starting the Mixing Process

Procedure

- ▶ If the “Mixing with recipe” or “Manual mixing” operating mode is selected: Press the [Start] button.
- ▶ If the “Remote mixing” operating mode is selected: Start the mixing process on the remote control unit (see the remote control unit instructions).
- ▷ The remaining time or the elapsed time for the mixing process or for a phase in a mixing process is displayed.

8.9.2 Pausing, Stopping, or Canceling the Mixing Process

Procedure

- ▶ If the mixing process in the “Mixing with recipe” or “Manual mixing” operating mode needs to be paused or stopped: Press the [Pause] or [Stop] button.
- ▶ If the mixing process in the “Remote mixing” operating mode needs to be paused: Pause the mixing process on the remote control unit (see the remote control unit instructions).
- ▶ If the mixing process has been stopped:
 - ▷ The [Stop mixing] display appears.
 - ▶ To confirm the process: Press the [YES] button.
 - ▷ The [Separate motor] display appears. The device must be decoupled from the Palletank® for Mixing.

8.10 Processing Messages

If faults occur during the mixing process when coupling or decoupling the device to | from the impeller and Palletank® for Mixing: A message appears in the status bar and the device gives off an acoustic signal.

Messages must be confirmed. All processes carried out on the device are interrupted until the messages are confirmed.

Requirements

An error message or warning message is displayed in the status bar.

Procedure

- ▶ Press the [Error] or [Warning] button.
- ▷ The error list appears.
- ▶ To switch between the entries in the error list: Press the [List up] or [List down] button.
- ▶ Check the entries in the error list:

- ▶ To call up the information for a message: Press the [Info] button.

- ▶ To call up the details for a message: Press the [Details] button.
- ▶ Resolve the fault, that has triggered the error message or warning message (see Chapter "10 Malfunctions", page 49).
- ▶ If the fault has been corrected: Confirm the message. To do so, press the [Confirm] button.
- ▷ The mixing process continues.

8.11 Viewing Recorded Messages

All recorded messages are saved. The in and out times of a message are marked by a time stamp. If the message memory is full: The oldest messages are deleted.

Procedure

- ▶ Call up the "System information" menu in the administration menu.
- ▶ Press the [Error history] button.
- ▷ The messages received are shown in a list:
 - Status K: In time of the message
 - Status KG: Out time of the message after confirmation by the user
- ▶ To switch between the entries in the error list: Press the [List up] or [List down] button.

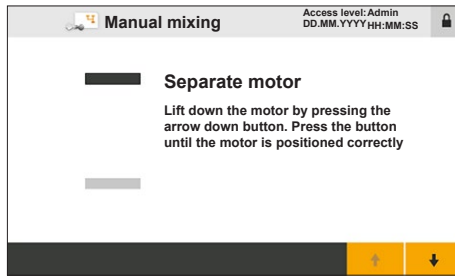
8.12 Ending the Mixing Process

8.12.1 Decoupling the Device from the Impeller and Palletank® for Mixing

To decouple the device from the Palletank® for mixing and the impeller: The motor, the alignment pins and the distance sensor must be lowered into the original position.

Requirements

- The mixing process is complete.
- The "Separate motor" display appears in the user guide.

Procedure

- ▶ Check the position of the motor in the “Separate motor” display. The positioning bars must be in the top position (1).
- ▶ Click on the [Lower] button and hold it down.
- ▶ The device is disconnected from the Palletank® for Mixing and the impeller in the mixing bag. During this process the distance sensor, the alignment pins, and the motor are lowered into the starting position and the device gives off acoustic signals.

8.12.2 Emptying and Removing the Mixing Bag

The mixing bag in the Palletank® for Mixing needs to be emptied and removed from the Palletank® for Mixing (see instructions for the Palletank® for Mixing).

8.12.3 Removing the Device from the Palletank® for Mixing

Requirements

The device is decoupled from the impeller and Palletank® for Mixing.

⚠ WARNING

Risk of injury when removing the device from the Palletank® for Mixing!

- ▶ Ensure that there are **no** people in the operating area of the device or the Palletank® for Mixing.

⚠ CAUTION

Risk of crushing due to moving parts!

When removing the device from the Palletank® for Mixing, the device’s guide rollers glide over the guide rails of the Palletank® for Mixing.

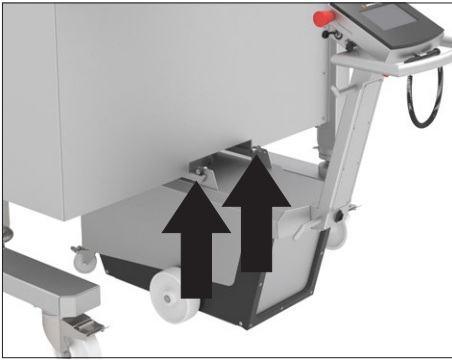
- ▶ Do **not** reach into the area between the guide rollers and guide rails of the Palletank® for Mixing.
- ▶ Ensure that there are **no** people in the operating area of the device who can reach between the guide rollers and guide rails.

⚠ CAUTION


Trip hazard from connection cable!

- ▶ While moving the device pay attention to where the connection cable is.
-

Procedure

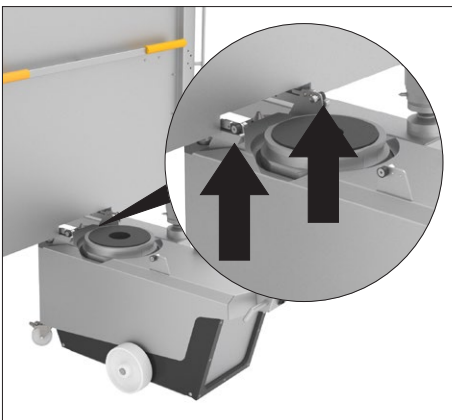- ▶ Slowly pull the device away from the Palletank® for mixing until the rear guide rollers are at the end of the guide rails of the Palletank® for mixing. The device must roll away over the 2 rolling elements on the guide rails of the Palletank® for Mixing.
- ▶ Pull the handle upwards slightly so that the rear guide rollers glide downwards from the guide rails of the Palletank® for Mixing.

- ▶ Continue to slowly pull the device away from the Palletank® for mixing until the front guide rollers are at the end of the guide rails of the Palletank® for mixing. The device must roll away over the 2 rolling elements on the guide rails of the Palletank® for Mixing.

- ▶ Press the handle downwards slightly so that the front guide rollers glide from the guide rails of the Palletank® for Mixing.
- ▶ If the device was used at a lower temperature than the temperature at the installation site where the device is intended to be used next: Acclimatize the device (see Chapter "6.2", Page31).
- ▶ **⚠ WARNING** Risk of injury from power supply when **not** in use! If the device is temporarily **not** in use: Switch off the device. Disconnect the power supply at the installation site and secure the power supply cable to the device's cable reel.

8.13 Ending the Process Data Acquisition

Requirements

The mixing process is complete.

Procedure

- ▶ Remove the USB stick from the device.
- ▶ Seal the connection with the cover cap.

9 Cleaning and Maintenance

9.1 Cleaning Device Surfaces

Requirements

The device is removed from the Palletank® for Mixing (see Chapter “8.12 Ending the Mixing Process”, page 45).

Procedure

- ▶ **⚠ CAUTION** Risk of injury due to uncontrolled rolling away of the device! Place it on a level surface and secure the brakes. Wear personal protective equipment.
- ▶ Turn the device off.
- ▶ Disconnect the power supply cable from the power supply at the installation site.
- ▶ If the device is connected to an external control system or other system: Remove the connection cable and close the connection with the closing cap.
- ▶ **NOTICE** Corrosion or damage to the device due to unsuitable cleaning agents!
 - ▶ Do **not** use corrosive, chloride-containing, and aggressive cleaning agents.
 - ▶ Do **not** use cleaning agents that contain abrasive ingredients, e.g. scouring agents, steel wool.
 - ▶ Do **not** use solvent-based cleaning agents.
 - ▶ Only use suitable cleaning agents (materials see Chapter “15.8 Materials”, page 59) and observe the product information for the cleaning agent used.
- ▶ **⚠ WARNING** Risk of injury due to hot surfaces! Allow the motor to cool down.
- ▶ Wipe the housing with a slightly damp cloth. Use a mild soapy solution or a suitable cleaning agent for more severe contamination.

9.2 Maintenance Schedule

Interval	Component	Action
12 months	Device	Have the function of the device checked by Sartorius Service.
12 months	Sensors	Have the device sensors calibrated by Sartorius Service.

9.3 Restarting the System

Procedure

- ▶ Commission the device (see Chapter “6 Getting Started”, page 30).

10 Malfunctions

10.1 Warning Messages

Warning Message	Malfunction (Info)	Cause	Solution	Chapter, Page
[Mixing warning: Mixing motor position]	[The measured position has changed more than 2 mm from the initial position]	During the mixing process, the distance sensor measures a signal that deviates from the output signal by more than 2 mm.	Check whether the impeller is moving correctly and freely (visual check in Palletank® for mixing). If required: Check the coupling of the device.	10.6, 52
[Mixing warning: Cart-in-place sensor]	[No cart-in-place sensor signal]	During the mixing process, the Palletank® recognition sensor loses the signal.	Check the coupling of the device.	10.6, 52
[Coupling warning: Impeller]	[No impeller detected in the mixing position]	The motor has reached the end position and cannot recognize the impeller.	Stop the mixing process. Decouple the device from the impeller and Palletank® for mixing. Check whether the impeller is correctly installed in the Palletank® for mixing (see instructions for the Palletank® for mixing). Couple the device with the impeller and Palletank® for mixing.	8.12.1, 45 8.6, 42
[Lifting system warning: Top end position reached]	[Lifting system top end position reached.]	The motor has reached the top end position and has not detected the alignment pins of the Palletank®.	Check that the distance sensor on the device or the alignment pin of the Palletank® is correctly adjusted (see Palletank® guide).	8.5, 40

10.2 Error Messages

Error Message	Malfunction (Info)	Cause	Solution	Chapter, Page
[Lifting system error: Motor timeout]	[Lifting motor does not reach position. (Sensor or motor are damaged)]	The motor cannot be correctly started-up, because the motor and other components are damaged.	Please contact Sartorius Service.	

Error Message	Malfunction (Info)	Cause	Solution	Chapter, Page
[Lifting system error: Sensor]	[Position sensors are damaged or wrongly calibrated]	The motor cannot be correctly started up, because the sensors are damaged or incorrectly calibrated.	Please contact Sartorius Service.	
[Mixing motor error: Temperature motor]	[Motor temperature exceeds limit (> 90°C)]	The temperature of the device's motor is too high, causing the mixing process to be interrupted and locked.	Ensure that the ambient temperature is within the acceptable range. If the error message reoccurs: Please contact Sartorius Service.	15.2, 57
[Mixing motor error: Temperature controller]	[Controller temperature exceeds limit (> 65°C)]	The temperature of the temperature regulator is too high, causing the mixing process to be interrupted and locked.	Ensure that the ambient temperature is within the acceptable range. If the error message reoccurs: Please contact Sartorius Service.	15.2, 57
[Mixing motor error: Mains voltage]	[Line voltage supply out of range or interrupted.]	The device is not being supplied with power or not sufficient power.	Check the power supply at the installation site.	15.3.1, 57
[Mixing motor error: Controller type]	[Motor not recognized.]	The connection to the motor is defective.	Please contact Sartorius Service.	
[Mixing motor error: Sensor cable]	[No sensor cable connected]	The connection to the motor is defective.	Please contact Sartorius Service.	
[Mixing motor error: Power cable]	[No power cable connected (or open bearing coil)]	The power supply cable is not connected or improperly connected.	Switch off the device. Check whether the power supply cable is properly connected. Switch on the device.	6.1.2, 30
[Mixing motor error: Drive current]	[Drive currents exceed limit (at least one phase > 11 A rms)]	The drive current is too high, causing the mixing process to be interrupted and locked.	Check the position of the mixing bag and impeller.	10.5, 52

Error Message	Malfunction (Info)	Cause	Solution	Chapter, Page
[Mixing motor error: Bearing current]	[Power consumption too high]	The storage current is too high, causing the mixing process to be interrupted or locked.	Check the position of the mixing bag and impeller.	10.5, 52
[Mixing motor error: Impeller unstable]	[Impeller is unstable]	The impeller in the mixing bag is unstable, causing the mixing process to be interrupted or locked.	Check the position of the mixing bag and impeller.	10.5, 52
[Mixing motor error: Drive coil]	[Open drive coil]	The Motor has a fault, causing the mixing process to be interrupted and locked.	Switch off the device. Check whether the power supply cable is properly connected. Switch on the device.	6.1.2, 30
[Mixing motor error: Power output]	[Controller power output short circuit or overtemperature]	The listed output is exceeded, causing the mixing process to be interrupted and locked.	Switch the device off and on again.	
[Mixing motor error: Speed lost]	[The rpm of the impeller is 0]	The mixing speed decreases to 0 during the mixing process and does not increase again within 10 seconds.	Stop the mixing process. Decouple the device from the impeller and Palletank® for mixing. Couple the device with the impeller and Palletank® for mixing.	8.12.1, 45 8.6, 42
[Mixing motor error: No impeller detected.]	[No impeller in the drive - either not present or too far away.]	The impeller is not recognized.	Check the position of the mixing bag and impeller.	10.5, 52

10.3 Faults when Operating the Device

Fault	Cause	Correction	Chapter, page
The device will not turn on.	The emergency shut-off switch is engaged.	Disengage the emergency shut-off switch. To do so, pull out the emergency shut-off switch until the green ring on the emergency shut-off switch is visible.	
The program menu or the administration menu are not displayed.	The necessary authorizations have not been assigned to the personal user profile.	Check the authorizations of the user profile.	13.1, 55

10.4 Leaked Liquids

Fault	Cause	Correction	Chapter, page
The device has come into contact with liquid because large amounts of liquids have leaked out of the Palletank® for Mixing.	The mixing bag is damaged and has fully or partially emptied. The liquid may have penetrated into the device via the fan.	Contact Sartorius Service.	

10.5 Checking the Position of the Mixing Bag and Impeller

Procedure

- ▶ Decouple the device from the impeller and Palletank® for Mixing (see Chapter "8.12.1", Page 45).
- ▶ Check that the bag and the impeller are correctly installed in the Palletank® for Mixing (see instructions for the Palletank® for Mixing).
 - ▶ If required: Adjust the position of the mixing bag.
- ▶ Insert the device under the Palletank® for Mixing and couple with the mixing bag's impeller (see Chapter "8.5", Page 40 and Chapter "8.6", Page 42).
 - ▶ If the warning message reoccurs: Contact Sartorius Service.

10.6 Checking the Coupling with the Mixing Bag and Impeller

Procedure

- ▶ Stop the mixing process.
- ▶ Decouple the device from the impeller and Palletank® for Mixing (see Chapter Chapter "8.12.1", Page 45).
- ▶ Wipe the Palletank® recognition sensor on the device with a damp cloth.
- ▶ Insert the device under the Palletank® for Mixing and couple with the mixing bag's impeller (see see Chapter "10.5", Page 52 and Chapter "8.6", Page 42).

11 Decommissioning

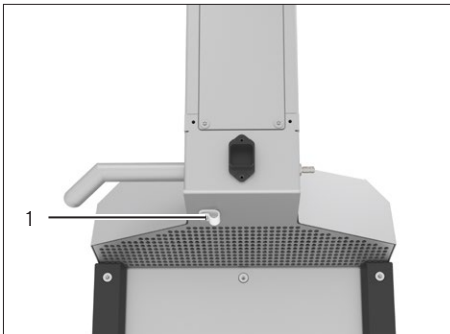
11.1 Decommissioning the Device

Requirements

The device is removed from the Palletank® (see Chapter “8.12 Ending the Mixing Process”, page 45).

Procedure

- ▶ Turn the device off.
- ▶ Disconnect the power supply cable from the power supply at the installation site.
- ▶ Disconnect the potential equalization cable from the power supply at the installation site.
- ▶ Unscrew the nut for the strain relief (1).
- ▶ Remove the power supply cable from the cable holder.
- ▶ Screw in the nut for the strain relief.
- ▶ Remove the potential equalization cable from the “Potential equalization” connection.
- ▶ If the device is connected to a remote control: Remove the connection cable and seal the connection with the locking cover cap.
- ▶ **⚠ CAUTION** Risk of injury due to uncontrolled rolling away of the device! Place it on a level surface and secure the brakes. Wear personal protective equipment.

12 Transport

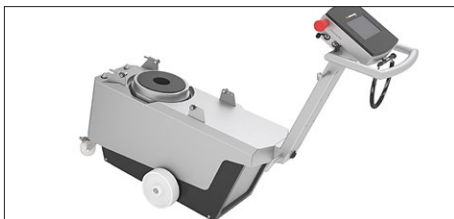
12.1 Installing the Transport Lock

If the device needs to be transported, e.g. if it needs to be shipped: The transport lock must be installed.


The transport lock consists of 2 transport lock screws that were fitted on the device on delivery.

Personnel: 2 persons, user


Procedure

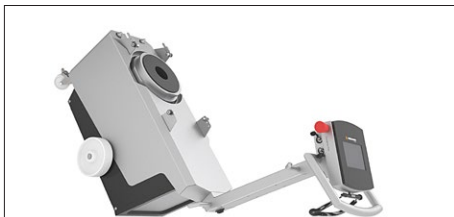- ▶ Hold the device firmly by the handle and push it down in a controlled manner.

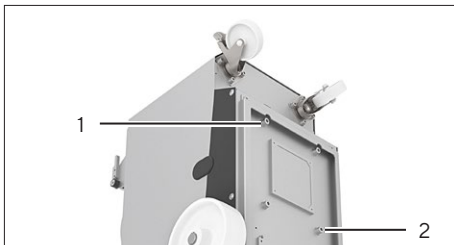- ▶ Push the device down until the handle is resting on the floor.
- ▶ Have a second person hold the device firmly by the handle, so the device remains in this position during installation.

- ▶ Unscrew screws (1) and (2), remove them, and save them for use with the device later.
- ▶ Screw the 2 transport lock screws into the device.
- ▶ While holding the device by the handle, tip it downwards once again in a controlled manner so that it stands on the castors.
- ▶ **⚠ CAUTION** Risk of injury due to uncontrolled rolling away of the device! Place it on a level surface and secure the brakes. Wear personal protective equipment.

13 Storage and Shipping

13.1 Storage

Procedure

- ▶ If the device is in operation:
 - ▶ Decommission the device.
 - ▶ Clean the device.
- ▶ Transport the device to the desired installation site (see Chapter “6.4 Moving the Device to the Installation Site”, page 32).
- ▶ Store the device according to the ambient conditions (see Chapter “15.2 Ambient Conditions”, page 57).
- ▶  **CAUTION** Risk of injury due to uncontrolled rolling away of the device! Place it on a level surface and secure the brakes. Wear personal protective equipment.

13.2 Returning Devices and Parts

Defective devices or parts can be returned to Sartorius Stedim Biotech. Returned devices must be clean, decontaminated, and properly packed.

Transport damage as well as measures for subsequent cleaning and disinfection of the device or parts by Sartorius Stedim Biotech shall be charged to sender.

WARNING

Risk of injury due to contaminated devices!

Devices contaminated with hazardous materials (nuclear, biological, or chemical NBC) will **not** be accepted for repair or disposal.

- ▶ Observe the information on decontamination (see Chapter “14.1”, Page 56).
-

Procedure

- ▶ Decommission the device.
- ▶ Install the transport lock.
- ▶ Contact Sartorius Service for instructions on how to return devices or parts (please refer to our website at www.sartorius.com for return instructions).
- ▶ Pack the device and its parts properly for return.

14 Disposal

14.1 Information on Decontamination

The device does **not** contain any hazardous materials that would necessitate special disposal actions.

The cultures and media (e.g., acids, bases) used during the processes are potentially hazardous materials that could cause biological or chemical hazards.

If the device has come into contact with hazardous substances: Steps must be taken to ensure proper decontamination and declaration. The operator is responsible for adhering to local government regulations on the proper declaration of transport and disposal and the proper disposal of the device.

14.2 Disposing of the Device and Parts

14.2.1 Information on Disposal

The device and the device accessories must be disposed of properly by disposal facilities.

The packaging is made of environmentally friendly materials that can be used as secondary raw materials.

14.2.2 Disposal

Requirements

The device has been decontaminated.

Procedure

- ▶ Dispose of the device. Follow the disposal instructions on our website (www.sartorius.com).
- ▶ Dispose of the packaging in accordance with local government regulations.

15 Technical Data

15.1 Dimensions and Weight

	Unit	Value
Dimensions (width x length x height)	mm	439 x 1242 x 1006
Weight	kg	100

15.2 Ambient Conditions

	Unit	Value
Installation Site		
Indoor use only		
Away from direct sunlight		
Not in close proximity to heating and air conditioning systems		
Away from radiation exposure from microwaves, UV lasers or X-radiation.		
Free from vibrations		
Maximum height above sea level	m	2000
Temperature	°C	+2 - +30
If temperature changes, i.e. due to changing locations, the device must be acclimatized		
Relative humidity, maximum		
At temperatures of up to 31 °C	%	80
Protection class according to IEC 60529: motor housing		IP 23
Protection class according to IEC 60529: electrical cabinet		IP 54
Pollution level according to IEC 61010-1		2

15.3 Electrical Data

15.3.1 Power Supply

	Unit	Value
Voltage	V _{AC}	100 - 240
Frequency	Hz	50/60
Performance (Current consumption), maximum	W	1200
Connection to the power supply		
Only use the original power supply cable supplied		
Length of the power supply cable	m	10

	Unit	Value
Minimum cross-sectional area of the connection leads for potential equalization	mm ²	2.5
Protection class according to IEC 60529		1

15.3.2 Electromagnetic Compatibility

	Unit	Value
Safety of electrical equipment, as per EN 61326-1 Electrical equipment for measurement, control, and laboratory use – EMC requirements – Part 1: General requirements (IEC 61326-1)		
Interference resistance: Suitable for use in industrial areas		
Transient emissions: Class		A
Class A are devices that are suitable for use in all other areas except residential areas and areas that are connected to a low voltage network that (also) supplies residential buildings.		

15.4 Suitability of Palletank® for Mixing and Mixing Bags

Only Sartorius Flexsafe® Pro Mixer Bag		
Filling volume of 50L to 1000L and a maximum of two connections on the front and underside		
Filling volume of 50L to 1000L and three or more connections on the front and underside		
Filling volume of 1500L to 3000L		
Only Sartorius Palletank® for Mixing		
Sizes 50L 100L 200L 400L 650L 1000L 1500L 2000L 2500L 3000L		
Approved for operation with Flexsafe® Pro Mixer Bag		

15.5 Mixing Process

	Unit	Value
Operating Modes		
Local operation (Manual, Recipe)		
Remote operation		
Controllable functions		
Starting, pausing, or stopping the mixing process		
Speed of the impeller	rpm	20 – 750

15.6 Remote Operation

	Unit	Value
Approved external control system		
Sartorius FlexAct® COM		
Communication connection		
Using a connection cable for remote operation		
Type of connection cable for remote operation: Shielded 8-pin plug, M12, A coded, screw assembly with M16 thread, 8 x 0.25 mm ²		
Input signals from remote control		
Target value for mixing speed at 0 – 750 rpm	V	0 – 10
Target value for mixing speed at 1 rpm	mV	13.3
Start mixing process: Relay status		Closed
Stop mixing process: Relay status		Open
Output signals of the device in remote operation		
Actual value for mixing speed at 0 – 750 rpm	mA	0 – 20
Actual value for mixing speed at 1 rpm	µA	21.3
Remote operation activated: Relay status		Closed
Remote operation deactivated (local operation): Relay status		Open
Warning message or error message present: Relay status		Closed
No warning message or error message on the device: Relay status		Open

15.7 Recipe Storage and Message Memory

	Unit	Value
Recipe storage		
Number of recipes, max		25
Number of phases in the recipe, max.		10
Message memory		
Number of saved messages (error messages, warning messages), max.		100

15.8 Materials

	Unit	Value
Device surfaces		
Stainless steel		316L
Glass bead blasted	Ra	1.2

	Unit	Value
Operating display: Foil coated with polyester foil		
Castors: Polyamide		

15.9 Acoustic Emission

	Unit	Value
Acoustic pressure level, max.	dB (A)	60

16 Accessories

This table contains an excerpt of the accessories that can be ordered. For information on other products, contact Sartorius Stedim Biotech.

Item	Quantity	Order number
Country-specific power supply cable for		
EU	1	FMA303410
US	1	FMA303411
Switzerland	1	FMA303412
Australia	1	FMA303413
GB	1	FMA303415
Connection cable for remote operation, length 3 m	1	FMA303416
USB stick	1	1000061414

17 Sartorius Service

Sartorius Service is always available to answer any queries regarding the device. Please visit the Sartorius website (www.sartorius.com) for information about the service addresses, services provided, or to contact a local representative.

When contacting Sartorius Service with questions about a system or in the event of malfunctions, be sure to have the device information close at hand e.g., serial number, hardware, firmware, and configuration. Consult the information on the manufacturer's ID label and in the "System Information | Software" menu.

18 Conformity Documents

The attached documents confirm compliance of the device with the directives or standards cited.

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SARTORIUS

EC / EU Declaration of Conformity

Manufacturer Sartorius Stedim Systems GmbH
34302 Guxhagen, Germany

declares under sole responsibility that the equipment

Device type Flexsafe® Pro Mixer

Type series Palletank for Mixing | Flexsafe® Pro Mixer Bag | Powder Bag Holder | Pro Mixer Drive Unit

Model **Palletank for Mixing:** FXC301930¹ ... FXC301949¹, FXC301951¹ ... FXC301956¹,
FXC301958¹ ... FXC301971¹
+ Flexsafe® Pro Mixer Bag: FMS500000 ... FMS500139
+ Powder Bag Holder: FXA304216; FXA304217; FXA304218
Pro Mixer Drive Unit: FMD300001

in the form as delivered fulfils all the relevant provisions of the following European Directives and meets the applicable requirements of the harmonized European Standards including any amendments valid at the time this declaration was signed listed below:

	EMC ²	RoHS	Machines
Directive	2014/30/EU	2011/65/EU	2006/42/EG 2006/42/EC
Standard(s)	EN 61326-1:2013	EN 50581:2012	EN ISO 12100:2010 EN 61010-1:2010 ³

The person authorised to compile the technical file:

Sartorius Stedim Systems GmbH
Product Compliance
34302 Guxhagen, Germany

Sartorius Stedim Systems GmbH
Guxhagen, 2021-02-18

Ariane Rindlisbacher
Manager Quality

Volker Otto
Head of Operations Guxhagen

- 1: For devices that have a winch or/and can be retrofitted with a winch
- 2: For devices with a weighing system or/and Pro Mixer Drive Unit
- 3: Applied standard, which however is not harmonized for machines

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Geräteart Flexsafe® Pro Mixer

Baureihe Palletank for Mixing | Flexsafe® Pro Mixer Bag | Powder Bag Holder | Pro Mixer Drive Unit

Modell
Palletank for Mixing: FXC301930¹ ... FXC301949¹, FXC301951¹ ... FXC301956¹,
FXC301958¹ ... FXC301971¹
+ Flexsafe® Pro Mixer Bag: FMS500000 ... FMS500139
+ Powder Bag Holder: FXA304216; FXA304217; FXA304218
Pro Mixer Drive Unit: FMD300001

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Type d'appareil Flexsafe® Pro Mixer

Série Palletank for Mixing | Flexsafe® Pro Mixer Bag | Powder Bag Holder | Pro Mixer Drive Unit

Modèle **Palletank for Mixing :** FXC301930¹ ... FXC301949¹, FXC301951¹ ... FXC301956¹,
FXC301958¹ ... FXC301971¹
+ Flexsafe® Pro Mixer Bag : FMS500000 ... FMS500139
+ Powder Bag Holder : FXA304216 ; FXA304217 ; FXA304218
Pro Mixer Drive Unit : FMD300001

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Tipo de aparato Flexsafe® Pro Mixer

Serie Palletank for Mixing | Flexsafe® Pro Mixer Bag | Powder Bag Holder | Pro Mixer Drive Unit

Modelo **Palletank for Mixing:** FXC301930¹ ... FXC301949¹, FXC301951¹ ... FXC301956¹, FXC301958¹ ... FXC301971¹
+ Flexsafe® Pro Mixer Bag: FMS500000 ... FMS500139
+ Powder Bag Holder: FXA304216; FXA304217; FXA304218
Pro Mixer Drive Unit: FMD300001

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Serie Pallettank for Mixing | Flexsafe® Pro Mixer Bag | Powder Bag Holder | Pro Mixer Drive Unit

Modello **Pallettank for Mixing:** FXC301930¹ ... FXC301949¹, FXC301951¹ ... FXC301956¹,
FXC301958¹ ... FXC301971¹
+ Flexsafe® Pro Mixer Bag: FMS500000 ... FMS500139
+ Powder Bag Holder: FXA304216; FXA304217; FXA304218
Pro Mixer Drive Unit: FMD300001

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Ref. Certif. No.

DE 3 - 31047

IEC SYSTEM FOR MUTUAL RECOGNITION OF TEST CERTIFICATES FOR ELECTRICAL EQUIPMENT (IECEE) CB SCHEME

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Name and address of the applicant

Sartorius Stedim Systems GmbH
Robert-Bosch-Str. 5 - 7
34302 Guxhagen
GERMANY

Name and address of the manufacturer

Sartorius Stedim Systems GmbH
Robert-Bosch-Str. 5 - 7, 34302 Guxhagen, GERMANY

Name and address of the factory

Baumüller Dravinja d.o.o.
Delavska Cesta 10, Celje, 3210 Slovenske Konjice, SLOVENIA

Ratings and principal characteristics

Rated voltage: 100-240 Vac
Rated frequency: 50-60 Hz
Rated power: 1200 W
Protection class: I

Trade mark (if any)

Sartorius Stedim

Model/type Ref.

Pro Mixer Drive Unit

A sample of the product was tested and found to be in conformity with

IEC 61010-1:2010
IEC 61010-2-051:2015

as shown in the Test Report Ref. No. which forms part of this certificate

028-713123890-000

This CB Test Certificate is issued by the National Certification Body
Ce Certificat d'essai OC est établi par l'Organisme **National de Certification**

CB 073361 0089 Rev. 00

Date, 2019-09-10

Page 1 of 1

(Ralph Fischer)

TÜV SÜD Product Service GmbH • Certification Body • Ridlerstraße 65 • 80339 Munich • Germany

Product Service

TÜV SÜD
ZERTIFIKAT ◆ CERTIFICATE ◆ 認證證書 ◆ CERTIFICADO ◆ CERTIFIKAT ◆ CERTIFICATE

A4 / 07.17

CERTIFICATE

No. U8 073361 0088 Rev. 00

Holder of Certificate: Sartorius Stedim Systems GmbH
Robert-Bosch-Str. 5 - 7
34302 Guxhagen
GERMANY

Production Facility(ies): 105452
Certification Mark:

Product: Laboratory equipment

Model(s): Pro Mixer Drive Unit

Parameters: Rated voltage: 100-240 Vac
Rated frequency: 50-60 Hz
Rated power: 1200 W
Protection class: I

Tested according to: UL 61010-1:2012/R:2016-04
CAN/CSA-C22.2 No. 61010-1:2012/U2:2016-04
UL 61010-2-051:2015
CAN/CSA-C22.2 No. 61010-2-051:2015

This product was voluntarily tested to the relevant safety requirements referenced on this certificate. It can be marked with the certification mark above. The mark must not be altered in any way. This product certification system operated by TÜV SÜD America Inc. most closely resembles system 3 as defined in ISO/IEC 17067. Certification is based on the TÜV SÜD "Testing and Certification Regulations". TÜV SÜD America Inc. is an OSHA recognized NRTL and a Standards Council of Canada accredited Certification body.

Test report no.: 028-713123890-000

Date, 2019-09-06

(Ralph Fischer)

Sartorius Stedim Systems GmbH
Robert-Bosch-Strasse 5-7
34302 Guxhagen, Germany

Phone: +49 5665 407 0
www.sartorius.com

The information and figures contained in these instructions correspond to the version date specified below.

Sartorius reserves the right to make changes to the technology, features, specifications and design of the equipment without notice.

Masculine or feminine forms are used to facilitate legibility in these instructions and always simultaneously denote the other gender as well.

Copyright notice:

This instruction manual, including all of its components, is protected by copyright. Any use beyond the limits of the copyright law is not permitted without our approval.

This applies in particular to reprinting, translation and editing irrespective of the type of media used.

© Sartorius Germany

Last updated:

03 | 2021